UNIVERSAL 320 UNIVERSAL 320 R

ⒹⒺ	Bedienungsanleitung	10
ⒺⒹ	Operating Instructions	35
ⒻⓇ	Mode d'emploi	59
ⒾⓉ	Istruzioni per l'uso	85

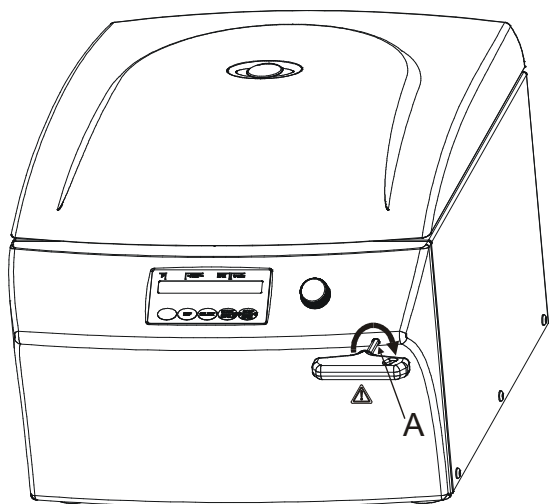
Fig. 1

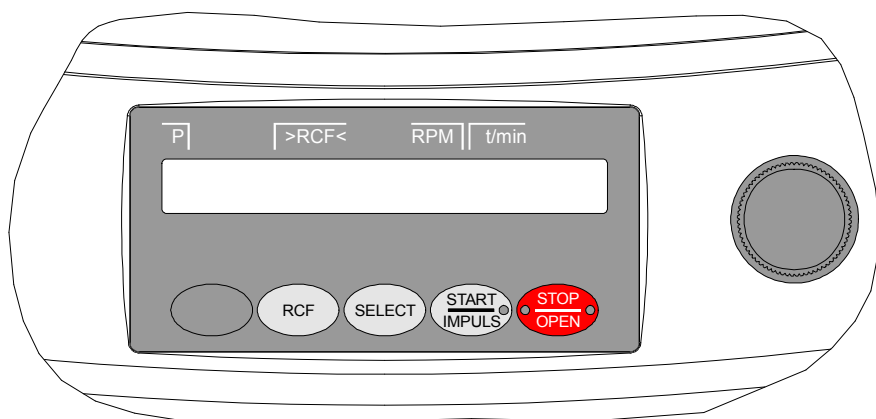
Fig. 2 UNIVERSAL 320

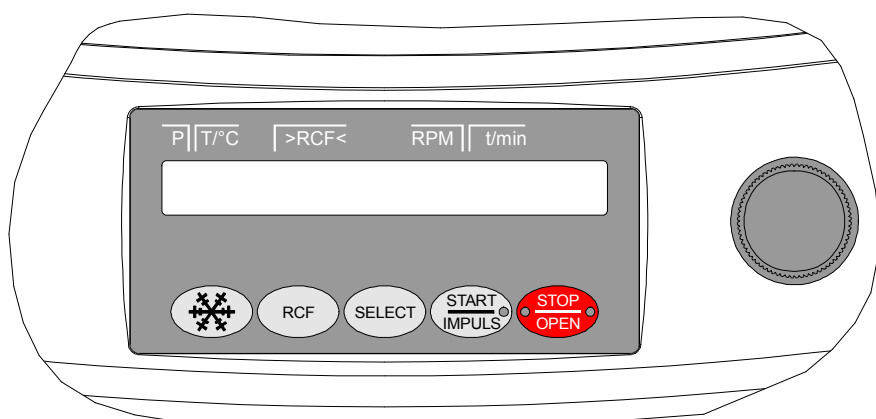
Fig. 3 UNIVERSAL 320 R

EG-Konformitätserklärung
EC Declaration of conformity
Déclaration de conformité CE
Dichiarazione di conformità CE

des Herstellers / of the manufacturer / du fabricant / del costruttore
Andreas Hettich GmbH & Co. KG • Föhrenstraße 12 • D-78532 Tuttlingen • Germany

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We hereby declare under our sole responsibility that the designated device and its accessories, which are listed in the technical documentation for this device and whose conformity has been assessed together with the device, conform to the Directive 98/79/EC on in vitro diagnostic medical devices.

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The conformity evaluation process was performed in accordance with appendix III of Directive 98/79/EC.

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The following additional European directives and ordinances have been applied:

- Machinery Directive 2006/42/EU
- EMC directive 2014/30/EU
- Low voltage directive 2014/35/EU
- RoHS II Directive 2011/65/EU (without involvement of a notified body)
- Ordinance (EC) No. 1907/2006 (REACH) (without involvement of a notified body)

Applied standards:

According to the list of applied standards, which is part of the product file.

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- **USA:** QSR, 21CFR 820 "CFR Title 21 - Food and Drugs: TITLE 21- FOOD AND DRUGS, CHAPTER I - FOOD AND DRUG ADMINISTRATION DEPARTMENT OF HEALTH AND HUMAN SERVICES, SUBCHAPTER H - MEDICAL DEVICES, Part 820 QUALITY SYSTEM REGULATIONS"
- **Kanada:** CMDR, SOR/98-282 "Medical Devices Regulations"

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Standards and regulations which apply to this device

The device is a high-end technical product. It is subject to extensive testing and certification procedures according to the following standards and regulations in their respectively valid version:

Electrical and mechanical safety for design and final testing:

Standard series: IEC 61010 (conform to standards of DIN EN 61010)

- IEC 61010-1 "Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements" (Pollution Degree 2, Excess-voltage category II)
- IEC 61010-2-010 "Safety requirements for electrical equipment for measurement, control and laboratory use - Part 2-010: Particular requirements for laboratory equipment for the heating of Materials" (only valid for centrifuges with heating)
- IEC 61010-2-011 "Safety requirements for electrical equipment for measurement, control and laboratory use - Part 2-011: Particular requirements for refrigerating equipment" (only valid for centrifuges with cooling)
- IEC 61010-2-020 "Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 2-020: Particular requirements for laboratory centrifuges"
- IEC 61010-2-101 "Safety requirements for electrical equipment for measurement, control and laboratory use - Part 2-101: Particular requirements for in vitro diagnostic (IVD) medical equipment"

Electromagnetic Compatibility:

- EN 61326-1 "Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements"

Risk management:

- DIN EN ISO 14971 "Application of risk management to medical devices"

Restriction of Hazardous Substances (RoHS II):

- EN 50581 "Technical documentation for assessing electric and electronic devices with regard to the restriction of hazardous substances"

European directives applied for conformity assessment procedures:

- In vitro diagnostic device directive 98/79/EG
EC conformity assessment procedure according to annex III "EC DECLARATION OF CONFORMITY" – self-declaration by the manufacturer
- Directive 2011/65/EU for the restriction of use of certain hazardous substances in electric and electronic devices. Carrying out the EC conformity assessment process is the sole responsibility of the manufacturer, without the involvement of a notified body.

Applied medical device regulations outside Europe:

- **USA:** QSR, 21CFR 820 "CFR Title 21 - Food and Drugs: TITLE 21- FOOD AND DRUGS, CHAPTER I - FOOD AND DRUG ADMINISTRATION DEPARTMENT OF HEALTH AND HUMAN SERVICES, SUBCHAPTER H - MEDICAL DEVICES, Part 820 QUALITY SYSTEM REGULATIONS"
- **Canada:** CMDR, SOR/98-282 "Medical Devices Regulations"

Certified quality management system according to

- ISO 9001 "Quality management systems – Requirements"
- ISO13485 "Medical devices - Quality management systems - Requirements for regulatory purposes"

Environmental management system according to

- ISO 14001 "Environmental management systems - Requirements with guidance for use"

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Restrictions relatives aux substances dangereuses (RoHS II):

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- **USA** : QSR, 21CFR 820 "CFR Title 21 - Food and Drugs : TITLE 21- FOOD AND DRUGS, CHAPTER I - FOOD AND DRUG ADMINISTRATION DEPARTMENT OF HEALTH AND HUMAN SERVICES, SUBCHAPTER H - MEDICAL DEVICES, Part 820 QUALITY SYSTEM REGULATIONS"
- **Canada**: CMDR, SOR/98-282 "Medical Devices Regulations"

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Contents

1	Use according to specification	37
2	Remaining risks	37
3	Technical specifications	37
4	Notes on safety	40
5	Transport and storage.....	41
5.1	Transport	41
5.2	Storage	41
6	Symbol meanings	41
7	Delivery checklist	42
8	Unpacking the centrifuge	42
9	Initial operation	43
10	Opening and closing the lid	43
10.1	Opening the lid	43
10.2	Closing the lid	43
11	Installation and removal of the rotor	44
11.1	Rotor with clamping nut.....	44
11.2	Rotor without clamping nut.....	44
11.2.1	Installing and removing the hub	44
11.2.2	Inserting and removing the rotor	44
12	Inserting and removing hangers into/from the rotor.....	45
13	Inserting and removing the adapters into/from the hangers	45
14	Inserting the adapter with positioning pin into the hanger	45
15	Loading the rotor	46
16	Handling of bio-safety systems.....	47
16.1	Opening and closing the lid of bio-safety systems	47
16.1.1	Lid with screw top and hole in turning handle	47
16.1.2	Lid with screw closure.....	48
16.1.3	Lid with clamp and tension lock	48
16.2	Storage of bio-safety systems	48
17	Control and display elements	48
17.1	Control knob.....	49
17.2	Control panel pushbuttons (keys)	49
17.3	Adjustment possibilities.....	49
18	Entering centrifugation parameter	50
19	Programming.....	50
19.1	Programme input/alteration.....	50
19.2	Programme recall.....	50
20	Centrifugation	50
20.1	Centrifugation with pre-set time	51
20.2	Continuous run.....	51
20.3	Short-term centrifugation.....	51
21	Emergency Stop.....	51
22	Acoustic Signal.....	51

23	Recall hours of operation	51
24	Cooling (only in centrifuges with cooling)	52
24.1	Standby-cooling	52
24.2	Pre-cooling the rotor	52
25	Relative centrifugal force (RCF)	52
26	Centrifugation of materials or mixtures of materials with a density higher than 1.2 kg/dm ³	53
27	Rotor recognition	53
28	Emergency release	53
29	Maintenance and servicing.....	54
29.1	Centrifuge (housing, lid and centrifuging chamber).....	54
29.1.1	Surface cleaning and care	54
29.1.2	Surface disinfection	54
29.1.3	Removal of radioactive contaminants	54
29.2	Rotors and Attachments	55
29.2.1	Cleaning and care.....	55
29.2.2	Disinfection	55
29.2.3	Removal of radioactive contaminants	55
29.2.4	Trunnions.....	56
29.2.5	Rotors and accessories with limited service lives	56
29.3	Autoclaving	56
29.4	Centrifuge containers.....	56
30	Faults	57
31	Returning Devices.....	58
32	Disposal	58
33	Anhang / Appen	111
33.1	Rotoren und Zubehör / Rotors and accessories.....	111

1 Use according to specification

This device is a medical product (laboratory centrifuge) within the context of the IVD Directive 98/79/EC.

The centrifuge was designed only for the separation of materials or mixtures with a density of no more than 1.2 kg/dm³. In particular, these include samples for preparation of in-vitro diagnostic purposes in human medicine.

The centrifuge is only meant for this purpose.

Another use or one which goes beyond this, is considered to be non-intended. The company Andreas Hettich GmbH & Co. KG is not liable for damage resulting from this.

Observing all information in the operating instructions and complying with the measures described therein is also a part of the intended use.

2 Remaining risks

The device is built according to the state-of-the-art and the recognized safety regulations. If used and handled improperly, there could be life-threatening danger to the user or third parties, or the device could be impaired or there could be other property damage. The device is only to be used for its intended purpose and only when it is in safe working condition.

Malfunctions which could affect safety must be corrected immediately.

3 Technical specifications

Manufacturer	Andreas Hettich GmbH & Co. KG D-78532 Tuttlingen				
Model	UNIVERSAL 320		UNIVERSAL 320 R		
Type	1401	1401-01	1406		1406-01
Mains voltage (± 10%)	200-240 V 1~	100-127 V 1~	200-240 V 1~	240 V 1~	115-127 V 1~
Mains frequency	50 – 60 Hz	50 – 60 Hz	50 Hz	60 Hz	60 Hz
Connected load	400 VA	400 VA	800 VA		950 VA
Current consumption	2.0 A	4.0 A	4.0 A		8.0 A
Cooling medium	----		R452A		
Max. capacity	4 x 200 ml				
Allowed density	1.2 kg/dm ³				
Speed (RPM)	16000				
Force (RCF)	24900				
Kinetic energy	9800 Nm				
Obligatory inspection (DGUV Regel 100 - 500)	no				
Ambient conditions (EN / IEC 61010-1) – Set-up site – Altitude – Ambient temperature – Humidity – Excess-voltage category (IEC 60364-4-443) – Pollution degree	Indoors only Up to 2000 m above sea level 2°C to 35°C 5°C to 35°C Maximum relative humidity 80% for temperatures up to 31°C, linearly decreasing to 50% relative humidity at 40°C. II 2				
Device protection class	I				
Not suitable for use in explosion-endangered areas.					
EMC – Emitted interference, Interference immunity	EN / IEC 61326-1, Class B	FCC Class B	EN / IEC 61326-1, Class B		FCC Class B
Noise level (dependent on rotor)	≤ 68 dB(A)		≤ 64 dB(A)		
Dimensions – Width – Depth – Height	401 mm 529 mm 346 mm		407 mm 698 mm 346 mm		
Weight	approx. 31 kg		approx. 52 kg		

Manufacturer	Andreas Hettich GmbH & Co. KG D-78532 Tuttlingen			
Model	UNIVERSAL 320 R			
Type	1406-50	1406-51	1406-70	1406-71
Mains voltage (± 10%)	200-240 V 1~	100-127 V 1~	200-240 V 1~	100-127 V 1~
Mains frequency	50 – 60 Hz	50 – 60 Hz	50 – 60 Hz	50 – 60 Hz
Connected load	400 VA	400 VA	400 VA	400 VA
Current consumption	2.0 A	4.0 A	2.0 A	4.0 A
Max. capacity	4 x 200 ml			
Allowed density	1.2 kg/dm ³			
Speed (RPM)	16000			
Force (RCF)	24900			
Kinetic energy	9800 Nm			
Obligatory inspection (DGUV Regel 100 - 500)	no			
Ambient conditions (EN / IEC 61010-1) – Set-up site – Altitude – Ambient temperature – Humidity – Excess-voltage category (IEC 60364-4-443) – Pollution degree	Indoors only Up to 2000 m above sea level 2°C to 35°C Maximum relative humidity 80% for temperatures up to 31°C, linearly decreasing to 50% relative humidity at 40°C. II 2			
Device protection class	I			
Not suitable for use in explosion-endangered areas.				
EMC – Emitted interference, Interference immunity	EN / IEC 61326-1, Class B	FCC Class B	EN / IEC 61326-1, Class B	FCC Class B
Noise level (dependent on rotor)	≤ 64 dB(A)			
Dimensions – Width – Depth – Height	401 mm 549 mm 346 mm		401 mm 550 mm 365 mm	
Weight	approx. 35 kg			

Manufacturer	Andreas Hettich GmbH & Co. KG D-78532 Tuttlingen		
Model	UNIVERSAL 320 R		
Type	1406-20		1406-21
Mains voltage (± 10%)	200-240 V 1~	240 V 1~	115-127 V 1~
Mains frequency	50 Hz	60 Hz	60 Hz
Connected load	800 VA		950VA
Current consumption	4.0 A		8.0 A
Cooling medium	R452A		
Max. capacity	4 x 200 ml		
Allowed density	1.2 kg/dm ³		
Speed (RPM)	16000		
Force (RCF)	24900		
Kinetic energy	9800 Nm		
Obligatory inspection (DGUV Regel 100 - 500)	no		
Ambient conditions (EN / IEC 61010-1) – Set-up site – Altitude – Ambient temperature – Humidity – Excess-voltage category (IEC 60364-4-443) – Pollution degree	Indoors only Up to 2000 m above sea level 5°C to 35°C Maximum relative humidity 80% for temperatures up to 31°C, linearly decreasing to 50% relative humidity at 40°C. II 2		
Device protection class	I		
Not suitable for use in explosion-endangered areas.			
EMC – Emitted interference, Interference immunity	EN / IEC 61326-1, Class B		FCC Class B
Noise level (dependent on rotor)	≤ 64 dB(A)		
Dimensions – Width – Depth – Height	407 mm 725 mm 365 mm		
Weight	approx. 52 kg		

4 Notes on safety

No claim of warranty will be considered by the manufacturer unless ALL instructions in this manual have been followed.

- The centrifuge should be installed on a good, stable base.
- Before using the centrifuge absolutely check the rotor for firm placement.
- When the centrifuge is running, according to EN / IEC 61010-2-020, no persons, dangerous substances or objects may be within the safety margin of 300 mm around the centrifuge.
- Rotors, suspensions and accessories that possess traces of corrosion or mechanical damage or if their term of use has expired may not be used any longer.
- The centrifuge may no longer be put into operation when the centrifuging chamber has safety-related damages.
- With swing-out rotors the trunnions must be regularly lubricated (Hettich Lubricating Grease No. 4051) in order to ensure consistent swinging out of the hangers.
- For centrifuges without temperature control, when the room temperature is increased and/or if the device is frequently used, the centrifuging chamber could be heated up. Therefore, it can't be ruled out that the sample material might be changed due to the temperature.

- Before the initial operation of your centrifuge you should read and pay attention to the operating instructions. Only personnel that has read and understood the operating instructions are allowed to operate the device.
- Along with the operating instructions and the legal regulations on accident prevention, you should also follow the recognised professional regulations for working in a safe and professional manner. These operating instructions should be read in conjunction with any other instructions concerning accident prevention and environmental protection based on the national regulations of the country where the device is to be used.
- This centrifuge is a state-of-the-art piece of equipment which is extremely safe to operate. However, it can lead to danger for users or others if used by untrained staff, in an inappropriate way or for a purpose other than that it was designed for.
- The centrifuge must not be moved or knocked during operation.
- In case of fault or emergency release, never touch the rotor before it has stopped turning.
- To avoid damage due to condensate, when changing from a cold to a warm room the centrifuge must either heat up for at least 3 hours in the warm room before being connected to the mains, or run hot for 30 minutes in the cold room.
- Only the rotors and accessories approved by the manufacturer for this device may be used (see chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories"). Before centrifuge vessels are used which are not listed in the chapter "Appendix, Rotors and accessories", the user must make sure they can be used by asking the manufacturer.
- The centrifuge rotor may only be loaded in accordance with the chapter "Loading the rotor".
- When centrifuging with maxim revolutions per minute the density of the materials or the material mixtures may not exceed 1.2 kg/dm³.
- The centrifuge may only be operated when the balance is within the bounds of acceptability.
- The centrifuge may not be operated in explosion-endangered areas.
- The centrifuge must not be used with:
 - inflammable or explosive materials
 - materials that react with one another producing a lot of energy.

- When centrifuging hazardous substances or mixtures, which are toxic, radioactive or contaminated with pathogenic microorganisms, suitable measures must be taken by the user.
Fundamentally, centrifuge containers with special screw closures must be used for hazardous substances. For materials of risk groups 3 and 4, in addition to sealable centrifuge containers, a bio-safety system must be used (see the "Laboratory Bio-safety Manual" from the World Health Organization).
In a bio-safety system, a bio-seal (sealing ring) prevents droplets and aerosols from escaping.
If the hanger of a bio-safety system is used without the lid, the sealing ring must be removed from the hanger to prevent damage to the sealing ring during the centrifugation run.
Damaged bio-safety systems are no longer microbiologically sealed.
If a bio-safety system is not used, a centrifuge is not microbiologically sealed for the purposes of the standard EN / IEC 61010-2-020.
When closing a bio-safety system, follow the instructions in the chapter "Handling of bio-safety systems".
For the available bio-safety systems, see the chapter "Appendix, Rotors and accessories". If in doubt, you can get the information you need from the manufacturer.
- The centrifuge must not be operated with highly corrosive substances which could impair the mechanical integrity of rotors, hangers and accessories.
- Repairs must only be carried out by personnel authorised to do so by the manufacturer.
- Only original spare parts and original accessories licensed by the Andreas Hettich GmbH & Co. KG company are allowed to be utilised.
- The following safety regulations apply:
EN / IEC 61010-1 and EN / IEC 61010-2-020 as well as their national deviations.
- The safe operation and reliability of the centrifuge can only be guaranteed if:
 - the centrifuge is operated in accordance with the operating instructions,
 - the electrical installation on the site where the centrifuge is installed conforms to the demands of EN / IEC stipulations.
- Meeting the country-specific requirements concerning occupational safety with regard to the use of laboratory centrifuges at the workplaces provided for this purpose by the user is the responsibility of the user.

5 Transport and storage

5.1 Transport

Before transporting the device, the transport securing device must be installed.

When the device and accessories are transported, the following ambient conditions must be complied with:

- Ambient temperature: -20°C to +60°C
- Relative humidity: 10% to 80%, non-condensing

5.2 Storage

The device and the accessories may only be stored in closed and dry rooms.

When the device and accessories are stored, the following ambient conditions must be complied with:

- Ambient temperature: -20°C to +60°C
- Relative humidity: 10% to 80%, non-condensing

6 Symbol meanings

Symbol on the device:
Attention, general hazard area.

Symbol on the device:
Observe operating instructions.
This symbol indicates that the user must observe the operating instructions provided.

Symbol in this document:

Attention, general hazard area.

This symbol refers to safety relevant warnings and indicates possibly dangerous situations.

The non-adherence to these warnings can lead to material damage and injury to personal.

Symbol on the device and in this document:

Beware of biohazard.

Symbol in this document:

This symbol refers to important circumstances.

Symbol on the device and in this document:

Symbol for the separate collection of electric and electronic devices according to the guideline 2002/96/EG (WEEE). The device belongs to Group 8 (medical devices).

Applies in the countries of the European Union, as well as in Norway and Switzerland.

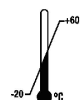
Symbol on the shipping carton label:

This way up.

Symbol on the shipping carton label:

The shipping packaging must be transported and handled within the indicated humidity range (10% - 80%).

Symbol on the shipping carton label:

The shipping packaging must be stored, transported and handled within the temperature range shown (-20°C - +60°C).

Symbol on the shipping carton label:

The shipping packaging must be kept away from rain and kept in a dry environment.

Symbol on the shipping carton label:

Fragile, handle with care.

7 Delivery checklist

- 1 Connecting cable
- 1 Hexagonal pin-type spanner
- 1 Lubricating grease for trunnions
- 1 Operating instructions
- 1 Notes on moving the equipment safely

The rotor(s) and associated accessories are included in the delivery in the quantity ordered.

8 Unpacking the centrifuge

- Lift the carton upward and remove the padding.

- 

Do not lift by the front panel.
Observe the weight of the centrifuge, refer to chapter "Technical specifications".

Lift the centrifuge on both sides with an appropriate number of helpers and place it on the laboratory table.

9 Initial operation

- Remove the transportation safety device from the bottom of the housing, see sheet "Transportation safety device"
- **Position the centrifuge in a stable and level manner in a suitable place. During set-up, the required safety margin of 300 mm around the centrifuge is to be kept according to EN / IEC 61010-2-020.**

When the centrifuge is running, according to EN / IEC 61010-2-020, no persons, dangerous substances or objects may be within the safety margin of 300 mm around the centrifuge.

- Ventilation openings may not be blocked.
A distance of 300 mm must be maintained from the ventilation slots and openings of the centrifuge.
- In the case of the centrifuge, type 1406-20, 1406-21, connect the nitrogen supply according to the enclosed instruction sheet AH1406-20XX.
In the case of the centrifuge, type 1406-50, 1406-51, connect the refrigerating/heating circulator according to the enclosed instruction sheet AH1406-50XX.
In the case of the centrifuge, type 1406-70, 1406-71, connect the refrigerating/heating circulator and the nitrogen supply according to the enclosed instruction sheet AH1406-70XX.

It is mandatory that the centrifuge be connected according to the enclosed instruction sheet.
Make sure you observe the enclosed instruction sheet.

- Check whether the mains voltage tallies with the statement on the type plate.
- Connect the centrifuge with the power cord to a standard mains socket. For connection ratings refer to Chapter "Technical specifications".
- Turn on the mains switch. Switch position "I".
The machine type and program version will be displayed and the LEDs light up. After 8 seconds, **OPEN** **OEFFNEN** is displayed and the left LED on the **STOP / OPEN** key blinks.
- Open the lid.
The last used centrifuge data will be displayed.

10 Opening and closing the lid

10.1 Opening the lid

The lid can only be opened when the centrifuge is switched on and the rotor is at rest. If it cannot be opened under these circumstances, see the section on "Emergency release".

- Press the button **OPEN / STOP**. The lid unlocks via the motor and the left LED in the pushbutton **OPEN / STOP** extinguishes.

10.2 Closing the lid

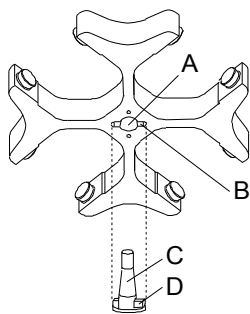
Do not put your fingers between lid and housing.
Do not bang the lid shut.

If the left LED in the **OPEN / STOP** button flashes, press the **OPEN / STOP** button so that the motor-driven lid lock goes into the basic position (opened).

- Place the lid and lightly press down the front edge of the lid. The locking action is effected by motor. The left LED in the button **OPEN / STOP** lights up.

11 Installation and removal of the rotor

11.1 Rotor with clamping nut

Rotor installation:

 Dirt particles between the motor shaft and rotor prevent the rotor from having a perfect seat and cause it to run unsteadily.

- Clean the motor shaft (C) and the bore of the rotor (A) and then apply a thin coat of grease to the motor shaft.
- Place the rotor vertically onto the motor shaft. The driver of the motor shaft (D) must be in the groove of the rotor (B). The groove alignment is marked on the rotor.
- Tighten the clamping nut of the rotor by turning it in the clockwise direction using the key included in delivery.
- Check the rotor to make sure it is seated firmly.

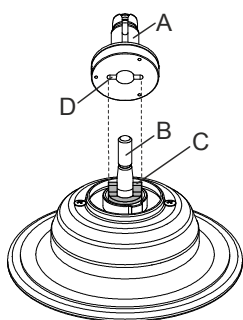
Removing the rotor:

- Loosen the clamping nut by turning it in the counterclockwise direction up to the lifting pressure point. After overcoming the lifting pressure point, the rotor is released from the cone of the motor shaft.
- Turn the clamping nut until the rotor can be lifted up from the motor shaft.

1.

11.2 Rotor without clamping nut

11.2.1 Installing and removing the hub

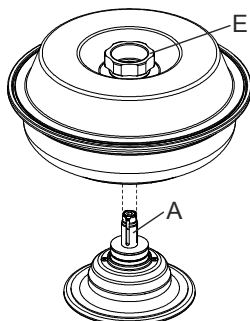
Installing the hub:

- Place the hub (A) vertically onto the motor shaft (B). The driver (C) on the motor shaft must be in the groove (D) of the hub.
- Tighten the clamping nut of the hub by turning it in the clockwise direction using the Allen key included in delivery.
- Check the hub to make sure it has a firm fit.

Removing the hub:

- Remove the rotor.
- Loosen the clamping nut of the hub by turning it counterclockwise up to the lifting pressure point. After overcoming the lifting pressure point, the hub is released from the cone of the motor shaft.
- Turn the clamping nut until the hub can be lifted up from the motor shaft.

11.2.2 Inserting and removing the rotor

Inserting the rotor:

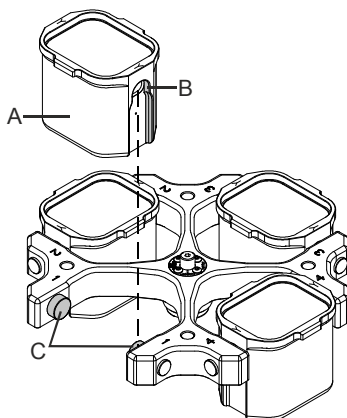
 Dirt particles between the hub and rotor prevent the rotor from having a perfect seat and cause it to run unsteadily. If necessary, clean the hub (A) on the motor shaft and the rotor hub before inserting the rotor.

- Lift up the rotor by the turning handle (E) of the lid, place it horizontally on the hub (A) and press it down as far as it can go.

Removing the rotor:

- Hold firmly onto the rotor by the turning handle (E) of the lid and lift it off of the hub (A).

12 Inserting and removing hangers into/from the rotor

For swing-out rotors, all rotor places must be occupied with the **same** hangers.

Certain hangers are marked with the number of the rotor place. These hangers may only be inserted in the corresponding rotor place.

Hangers which are labelled with a set number, e.g. S001/4, may only be used in a set.

Inserting hangers in the rotor:

- Check the rotor to make sure it is seated firmly.
- Grease the lifting lug (C) (Hettich lubricating grease, no. 4051).
- Insert hangers (A) into the rotor. While doing so, make sure that the lifting lugs (C) are in the grooves (B) of the hangers.
- Push the hangers downward as far as they can go.

Removing hangers from the rotor:

- Pull the hangers (A) straight upward and out of the rotor.

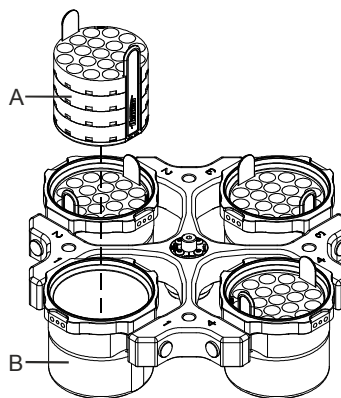
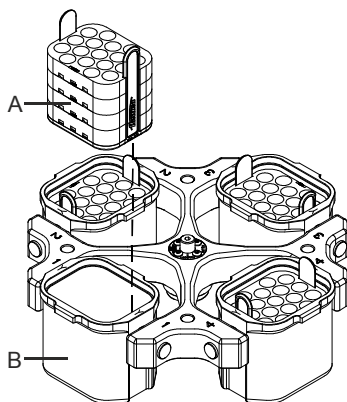
13 Inserting and removing the adapters into/from the hangers

Inserting the adapters into the hangers:

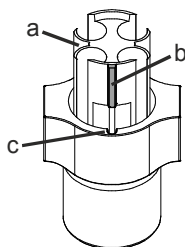
- Insert the adapters (A) horizontally into the hangers (B).

Remove the adapters from the hangers:

- Remove the adapters (A) upward and out of the hangers (B).

14 Inserting the adapter with positioning pin into the hanger

- Insert the adapter (a) into the hanger such that the positioning pin (b) is in the groove (c) of the hanger.

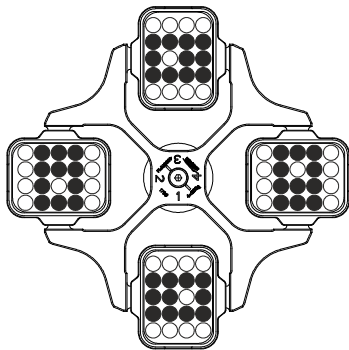
If the adapter (a) is not correctly inserted, the centrifuge containers could be damaged during the centrifugation run.

15 Loading the rotor

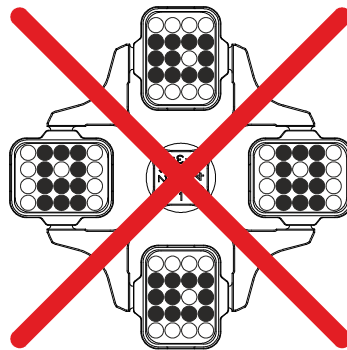
Standard centrifuge containers of glass will not stand RCF values exceeding 4000 (DIN 58970, pg. 2).

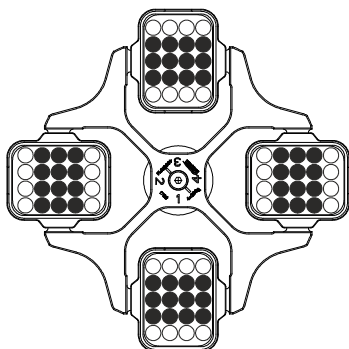
- Check the rotor for firm seating.
- With swing-out rotors all rotor positions must be lined with **identical** hangers. Certain hangers are marked with the number of the rotor position. These hangers may only be used in the respective rotor position. Hangers that are marked with a set number (e.g. S001/4) may only be used in the set.
- The rotors and hangers may only be loaded symmetrically. The centrifuge containers have to be distributed evenly on all rotor positions. For authorised combinations see Chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories".
In the case of angle rotors all possible rotor positions must be loaded, see chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories".

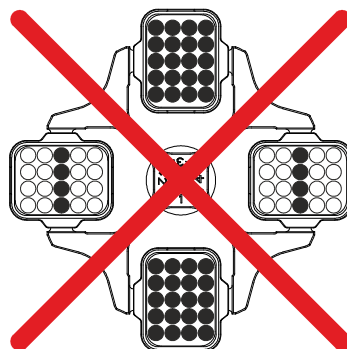
Rotor is loaded symmetrically.

Not permitted!
Rotor is loaded asymmetrically.

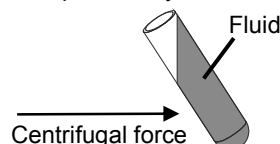
Rotor is evenly loaded

Not permitted!
Rotor is not evenly loaded

- On certain hangers, the weight of the maximum load or the weight of the maximum load and the maximum weight of the completely loaded hanger is specified. This weight may not be exceeded. In case of exception, see chapter "Centrifugation of materials or mixtures of materials with a density higher than 1.2 kg/dm³". The weight specification of the maximum load includes the total weight of the adapter, centrifuge container and contents.
- In containers with rubber inserts, the same number of rubber inserts must always be among the centrifuge containers.
- The centrifuge containers may only be filled outside of the centrifuge.

- The maximum filling quantity for the centrifuge containers specified by the manufacturer must not be exceeded.
In the case of angle rotors, the centrifuging vessels may only be filled so far that no fluid can be expelled from them while the centrifuge is running.

- When loading the angle rotors, no liquid may get into the angle rotors or centrifuging chamber.
- When loading the hangers and swing-out rotors and when swivelling out the hangers while the centrifuge is running, no liquid may enter the hangers or the centrifuging chamber.
- In order to maintain the weight differences within the centrifuge container as marginal as possible, a consistent fill level in the containers is to be heeded.

16 Handling of bio-safety systems

Bio-safety systems may only be used when they are dry.

Each time before using the bio-safety system, all parts of the bio-safety system must be visually inspected for damage. In addition, the sealing ring(s) of the bio-safety system must be checked to make sure they are in the correct installation position.

Damaged parts of the bio-safety system must be exchanged immediately.

Damaged bio-safety systems are no longer microbiologically sealed.

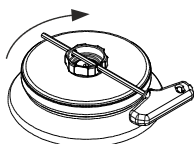
To ensure leak-tightness, the lid of the bio-safety system must be correctly sealed.

For the available bio-safety systems, see the chapter "Appendix, Rotors and accessories". If in doubt, you can get the information you need from the manufacturer.

16.1 Opening and closing the lid of bio-safety systems

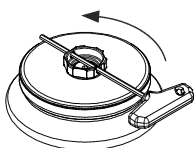
16.1.1 Lid with screw top and hole in turning handle

Closing the lid:

- Place the lid centred on the rotor.
- Stick the supplied key through the hole in the turning handle and close the lid firmly by turning the key in a clockwise direction.

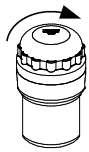
Opening the lid:

- Stick the supplied key through the hole in the turning handle and turn counter-clockwise until the lid can be removed from the rotor.

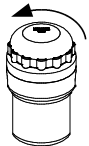
16.1.2 Lid with screw closure

Closing the lid:

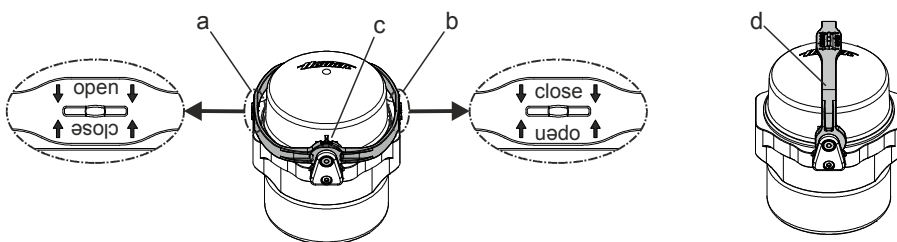
- Place the lid on the hanger.
- Tightly seal the lid by hand by turning it clockwise.

Opening the lid:

- Turn the lid by hand counterclockwise until the lid can be removed from the hanger.

16.1.3 Lid with clamp and tension lock

Closing the lid:

- Swivel the clamp in the "↓ open ↓" (a) position. The arrows of the "↓ open ↓" label must be pointing downwards so that the text "open" is legible.
- Place the lid on the hanger such that the two pins of the lid are in the two openings of the clamp (c).
- Swivel the clamp to the opposite side of the hanger to the "↓ close ↓" (b) position. The arrows of the "↓ close ↓" must be pointing downwards so that the text "close" is legible.

The clamp must lie on the hanger so that it can swing out during the centrifugation run.

- If the clamp is swivelled into position (d), the bio-safety system can be carried by the clamp.

The leak-tightness of the bio-safety system is also ensured in the carry position.

During transportation, do not swing the bio-safety system back and forth, since otherwise, leak-tightness can no longer be guaranteed.

- For easier handling when inserting and removing the hanger, swivel the clamp into position (d) and hold the hanger on the clamp.

Opening the lid:

- Swivel the clamp in the "↓ open ↓" (a) position. The arrows of the "↓ open ↓" label must be pointing downwards so that the text "open" is legible.
- Remove the lid from the hanger.

16.2 Storage of bio-safety systems

- To prevent damage to the sealing rings during storage, bio-safety systems may only be stored with the lid open.

1. Control and display elements

See figure on page 2.

Fig. 2, Fig. 3: Display and control panel

17.1 Control knob

For setting the individual parameters.
Turning anticlockwise reduces the value. Turning clockwise increases the value.

1.2 Control panel pushbuttons (keys)

- Selection control key for selection of specific parameter.
The subsequent parameter is selected by every further keystroke.

- Start centrifugation run. The LED in the button lights up during the centrifugation run as long as the rotor is turning.
- Short-term centrifugation.
The centrifugation run is effected as long as the button is held down. The LED in the button lights up during the centrifugation run as long as the rotor is turning.
- Store inputs and changes.

- End centrifugation run.
The rotor runs down with a pre-selected brake step. The right-hand LED in the button lights up until the rotor is stationary. Once the rotor is stationary the left-hand LED flashes in the button. Pressing the button twice triggers the EMERGENCY STOP.
- Unlock the lid.
The left-hand LED in the button goes out.
- Leave the parameter input.

- Switch between RPM and RCF display.
RCF values are displayed in > <.

- Start pre-cooling.
The pre-cooling speed is settable. It is pre-adjusted to 2800 RPM.

1.3 Adjustment possibilities

PROG RCL Program position of the called-up program.

t/min Running time. Settable from 0 - 99 min, in 1 min increments.

t/sec Running time. Settable from 0 - 59 s, in 1 second increments.

Continuous run "∞". Set parameter **t/min** and **t/sec** to zero.

RPM Revolutions per minute. A numerical value from 500 RPM up to the maximum speed of the rotor can be set. Maximum speed of the rotor, see Chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories". Settable in increments of 10.

RAD/mm Centrifugation radius. Input in mm. For centrifugation radius see Chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories". The input of the radius is only possible if the RCF display (> RCF <) is selected.

RCF Relative centrifugal force. A numerical value can be set, which gives a speed between 500 RPM and the maximum speed of the rotor. Adjustable up to 100 in intervals of 1, and from 100 in intervals of 10. The RCF value is automatically rounded up or rounded down with regard to the RPM interval. The input of the RCF is only possible if the RCF display (> RCF <) is selected.

Starting steps 1 - 9. Step 9 = shortest starting time, Step 1 = longest starting time.

Brake steps 0 - 9. Step 9 = shortest run-down time, Step 1 = long run-down time, Step 0 = longest run-down time (brakeless run-down).

T/°C Temperature Set Point (only in centrifuges with cooling). Adjustable from -20°C to +40°C, in 1°C intervals. The lowest obtainable temperature depends on the rotor (see Chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories").

PROG STO Program position on which the program is stored. 9 programs can be stored (program positions 1 - 2 - 3 - ... 9). The programme position # serves as temporary storage for altered adjustments.

1.

1. Entering centrifugation parameter

If no key is pressed for 8 seconds long after the selection or during the input of parameters, the previous values will be shown in the display. The input of parameter then has to be executed again.

- Select the RPM or RCF display with the key **[RCF]**. RCF values are displayed in $> <$.
- Select the desired parameters using the button **[SELECT]** and set using the knob $\odot$.
In order to set continuous operation, the parameters **t/min** and **t/sec** must be set to zero with the $\odot$ knob.
Continual running is represented in the display by the following symbol, " ∞ ".
- After input of all parameters, press the key **[START/IMPULS]** in order to store the adjustments on the programme position **#**. As confirmation, ***** ok ***** will be displayed for a short period.

The data on the programme position **#** will be overwritten with every input of parameters and pressing of the key **[START/IMPULS]**.

2. Programming

2.1 Programme input/alteration

If no key is pressed for 8 seconds long after the selection or during the input of parameters, the previous values will be shown in the display. The input of parameter then has to be executed again.

- Select the RPM or RCF display with the key **[RCF]**. RCF values are displayed in $> <$.
- Select the desired parameters using the button **[SELECT]** and set using the knob $\odot$.
In order to set continuous operation, the parameters **t/min** and **t/sec** must be set to zero with the $\odot$ knob.
Continual running is represented in the display by the following symbol, " ∞ ".
- The parameter **PROG STO** can be selected using the button **[SELECT]** and the desired program position set using the knob $\odot$.
- Press the button **[START/IMPULS]** in order to store the setting on the desired program position. ***** ok ***** is displayed briefly as confirmation.
If the key **[START/IMPULS]** is pressed without the parameter **PROG STO** being activated, the settings are always stored in the program place **#**.

The previous data in the program position is overwritten during saving.

2.2 Programme recall

- Select the parameter **PROG RCL** using the button **[SELECT]** and set the desired program position using the knob $\odot$.
- Press the button **[START/IMPULS]**. The centrifugation data of the selected program position is displayed.
- The parameters can be checked by pressing the button **[SELECT]**.
To leave the parameter display press the button **[OPEN/STOP]** or press no button for a period of 8 seconds.

3. Centrifugation

When the centrifuge is running, according to EN / IEC 61010-2-020, no persons, dangerous substances or objects may be within the safety margin of 300 mm around the centrifuge.

If the permissible weight difference within the rotor loading has been exceeded, the drive shuts down during the start-up, the unbalance display lights up, and **IMBALANCE** is displayed.

A centrifugation run can be stopped at any time by pushing the key **[OPEN/STOP]**.

All parameters can be selected and altered during the centrifugation run (see Chapter "Entering centrifugation parameter").

You can switch-over at any time between the RPM and RCF display with the key **[RCF]**. The input of the centrifugation radius is necessary if you are working with the RCF display.

If **OPEN** **OEFFNEN** is displayed, a further operation of the centrifuge is only possible after opening the lid once.

If **R xx n-max xxxxx** is displayed, then no centrifugation run has taken place as the rotor was changed beforehand, refer to Chapter "Rotor Identification".

- Turn on the mains switch. Switch position **I**.
- Load the rotor and close the centrifuge lid.

3.1 Centrifugation with pre-set time

- Adjusting time or recall a programme with pre-set time (see Chapter "Programming").
- Press the key **START/IMPULS**. The LED in the button **START/IMPULS** lights up for as long as the rotor turns.
- After expiration of the time or with truncation of the centrifugation run by pushing the key **OPEN/STOP**, the run-down is effected with the selected brake step. The brake step is displayed.

During the centrifugation run, the rotor speed or the RCF value, the temperature in the centrifuging chamber (only for centrifuge with cooling), and the remaining time are displayed.

3.2 Continuous run

- Adjusting the symbol ∞ or recall a continuous run programme (see Chapter "Programming").
- Press the key **START/IMPULS**. The LED in the button **START/IMPULS** lights up for as long as the rotor turns. The time metering begins at 00:00.
- Press the key **OPEN/STOP** in order to stop the centrifugation run. The run-down is effected with the selected brake step. The brake step is displayed.

During the centrifugation run, the rotor speed or the RCF value, the temperature in the centrifuging chamber (only for centrifuge with cooling), and the elapsed time are displayed.

3.3 Short-term centrifugation

- Hold down the key **START/IMPULS**. The LED in the button **START/IMPULS** lights up for as long as the rotor turns. The time metering begins at 00:00.
- Let go of the key **START/IMPULS** again in order to stop the centrifugation run. The run-down is effected with the selected brake step. The brake step is displayed.

During the centrifugation run, the rotor speed or the RCF value, the temperature in the centrifuging chamber (only for centrifuge with cooling), and the elapsed time are displayed.

4. Emergency Stop

- Press the key **OPEN/STOP** twice.

With Emergency Stop the run-down is effected with brake step 9 (shortest run-down time). Brake step 9 is displayed. If brake step 0 was pre-selected, the run-down time is technically longer than with brake step 9.

5. Acoustic Signal

The acoustic signal sounds:

- Upon the appearance of a disturbance in 2 second intervals.
- After completion of a centrifugation run and rotor standstill in 30 second intervals.

The acoustic signal is stopped by opening the lid or pressing any key.

The signal after completion of the centrifugation run can be activated or deactivated in the following manner, if the rotor is at standstill:

- Hold down the key **SELECT** for 8 seconds.
After 8 seconds, **SOUND / BELL** appears in the display.
- Set using the knob **OFF** or **ON**.
- Press the key **START/IMPULS** in order to store the setting.
As confirmation, ***** ok ***** will be displayed for a short period.

6. Recall hours of operation

Recall hours of operation is only possible during rotor standstill.

- Hold down the key **SELECT** for 8 seconds.
After 8 seconds, **SOUND / BELL** appears in the display.
- Press the key **SELECT** once again.
The centrifuge's hours of operation (**CONTROL:**) are displayed.
- Press the key **OPEN/STOP** to exit the hours of operation recall.

2.

24 Cooling (only in centrifuges with cooling)

The temperature set-point can be adjusted from -20°C to +40°C. The lowest obtainable temperature is dependent on the rotor (see Chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories").

24.1 Standby-cooling

With rotor standstill and closed lid the centrifugal chamber is cooled to the pre-selected temperature. The temperature set-point is shown in the display.

Standby cooling will be subject to a timed delay after a centrifuge run and the display will show **OPEN** **OEFFNEN**. The delay time can be pre-set in 1 minute steps from 1 to 5 minutes. It is pre-set to 1 minute.

With the rotor standing still and the cover open the delay time can be set as follows:

- Hold down the key  for 8 seconds.
After 8 seconds, **t/min = X** appears in the display.
- Use the rotary button  to set the delay time.
- Press the key  in order to store the setting.
As confirmation, ***** ok ***** will be displayed for a short period.

To leave the delay time display press the key  or do not press any key for a period of 8 seconds.

24.2 Pre-cooling the rotor

- Press the key . The LED in the button  lights up for as long as the rotor turns.
- Press the button  to end the pre-cooling. The run-down is effected with the selected brake step. The brake step is displayed.

During the centrifugation run, the rotor speed or the RCF value, the temperature in the centrifuging chamber, and the elapsed time are displayed.

The pre-cooling speed can be adjusted in decadic steps from 500 RPM to the max RPM of the rotor. It is pre-adjusted to 2800 RPM.

When the rotor is stationary and the lid open the pre-cooling speed can be set in the following manner:

- Hold down the key  for 8 seconds.
After 8 seconds, **t/min = X** appears in the display.
- Press the key  once again.
The set pre-cooling RPM - **RPM = XXXX** will be displayed.
- Set the desired pre-cooling speed using the knob .
- Press the key  in order to store the setting.
As confirmation, ***** ok ***** will be displayed for a short period.

To leave the pre-cooling RPM display press the key  or do not press any key for a period of 8 seconds.

25 Relative centrifugal force (RCF)

The relative centrifugal force (RCF) is given as a multiple of the acceleration of gravity (g). It is a unit-free value and serves to compare the separation and sedimentation performance.

These values are calculated using the formula below:

$$RCF = \left(\frac{RPM}{1000} \right)^2 \times r \times 1,118 \quad \Rightarrow \quad RPM = \sqrt{\frac{RCF}{r \times 1,118}} \times 1000$$

RCF = relative centrifugal force

RPM = rotational speed (revolutions per minute)

r = centrifugal radius in mm = distance from the centre of the turning axis to the bottom of the centrifuge container. For more on the centrifugal radius see the chapter "Anhang/Appendix, Rotoren und Zubehör/Rotors and accessories".

The relative centrifugal force (RCF) stands in relation to the revolutions per minute and the centrifugal radius.

26 Centrifugation of materials or mixtures of materials with a density higher than 1.2 kg/dm³

When centrifuging with maxim revolutions per minute the density of the materials or the material mixtures may not exceed 1.2 kg/dm³.

The speed must be reduced for materials or mixtures of materials with a higher density.

The permissible speed can be calculated using the following formula:

$$\text{Reduced speed (n}_{\text{red}}) = \sqrt{\frac{1.2}{\text{Greater density [kg/dm}^3]}} \times \text{maximum speed [RPM]}$$

e.g.: maximum speed RPM 4000, density 1.6 kg/dm³

$$n_{\text{red}} = \sqrt{\frac{1.2 \text{ kg/dm}^3}{1.6 \text{ kg/dm}^3}} \times 4000 \text{ RPM} = 3464 \text{ RPM}$$

In the exceptional case that the maximum loading indicated on the hanger is exceeded, the speed must also be reduced.

The permissible speed can be calculated using the following formula:

$$\text{Reduced speed (n}_{\text{red}}) = \sqrt{\frac{\text{maximum load [g]}}{\text{actual load [g]}}} \times \text{maximum speed [RPM]}$$

e.g.: maximum speed RPM 4000, maximum load 300 g, actual load 350 g

$$n_{\text{red}} = \sqrt{\frac{300 \text{ g}}{350 \text{ g}}} \times 4000 \text{ RPM} = 3703 \text{ RPM}$$

If in doubt you should obtain clarification from the manufacturer.

27 Rotor recognition

Rotor recognition is carried out after each start of the centrifugation run.

If the rotor has been changed, the centrifugation run is canceled after rotor recognition. The rotor code (R xx) and the maximum speed (n-max=xxxxx) of the rotor are displayed.

A further operation of the centrifuge is only possible after a single opening of the lid.

If the maximum speed of the rotor being used is less than the set speed, the speed is limited to the rotor's maximum speed.

1. Emergency release

During a power failure the lid cannot be unlocked by motor. An emergency release has to be executed by hand.

For emergency release disconnect the centrifuge from the mains.
Open the lid only during rotor standstill.

See figure on page 2.

- Switch off the mains switch (switch position "0").
- Look through the window in the lid to be sure that the rotor has come to a standstill.
- Insert the hexagonal wrench key into the bore hole (Fig. 1, A) and carefully rotate by half a turn in clockwise direction until the lid can be opened.
- Pull the hexagon socket head wrench out of the drilling again.
- Re-insert the plug.
- If the left LED in the **OPEN / STOP** button flashes after the centrifuge is switched on again, press the **OPEN / STOP** button so that the motor-driven lid lock goes into the basic position (opened) again.

29 Maintenance and servicing

The device can be contaminated.

Pull the mains plug before cleaning.

Before any other cleaning or decontamination process other than that recommended by the manufacturer is applied, the user has to check with the manufacturer that the planned process does not damage the device.

- Centrifuges, rotors and accessories must not be cleaned in rinsing machines.
- They may only be cleaned by hand and disinfected with liquids.
- The water temperature must be between 20 – 25°C.
- Only detergents/disinfectants may be used which:
 - have a pH between 5 - 8
 - do not contain caustic alkalis, peroxides, chlorine compounds, acids and alkaline solutions
- In order to prevent appearances of corrosion through cleaning agents or disinfectants, the application guide from the manufacturer of the cleaning agent or disinfectant are absolutely to be heeded.

29.1 Centrifuge (housing, lid and centrifuging chamber)

29.1.1 Surface cleaning and care

- Clean the centrifuge housing and the centrifuging chamber regularly, using soap or a mild detergent and a damp cloth if required. For one thing, this services purposes of hygiene, and it also prevents corrosion through adhering impurities.
- Ingredients of suitable detergents: soap, anionic tensides, non-ionic tensides.
- After using detergents, remove the detergent residue by wiping with a damp cloth.
- The surfaces must be dried immediately after cleaning.
- In the event of condensation water formation, dry the centrifugal chamber by wiping out with an absorbent cloth.
- Lightly rub the rubber seal of the centrifuge chamber with talcum powder or a rubber care product after each cleaning.
- The centrifuging chamber is to be checked for damage once a year.

If damage is found which is relevant to safety, the centrifuge may no longer be put into operation. In this case, notify Customer Service.

29.1.2 Surface disinfection

- If infectious materials penetrates into the centrifugal chamber this is to be disinfected immediately.
- Ingredients of suitable disinfectants: ethanol, n-propanol, ethyl hexanol, anionic tensides, corrosion inhibitors.
- After using disinfectants, remove the disinfectant residue by wiping with a damp cloth.
- The surfaces must be dried immediately after disinfecting.

29.1.3 Removal of radioactive contaminants

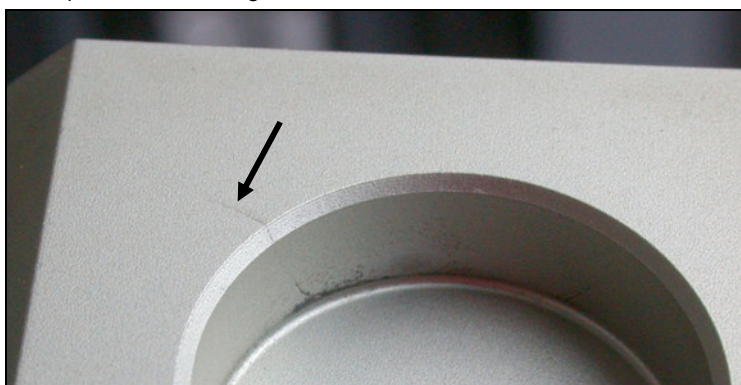
- The agent must be specifically labelled as being an agent for removing radioactive contaminants.
- Ingredients of suitable agents for removing radioactive contaminants: anionic tensides, non-ionic tensides, polyhydrated ethanol.
- After removing the radioactive contaminants, remove the agent residue by wiping with a damp cloth.
- The surfaces must be dried directly after removing the radioactive contaminants.

29.2 Rotors and Attachments

29.2.1 Cleaning and care

- In order to avoid corrosion and changes in materials, the rotors and accessories have to be cleaned regularly with soap or with a mild cleaning agent and a moist cloth. Cleaning is recommended at least once a week. Contaminants must be removed immediately.
- Ingredients of suitable detergents: soap, anionic tensides, non-ionic tensides.
- After using detergents, remove detergent residue by rinsing with water (only outside of the centrifuge) or wipe off with a damp cloth.
- The rotors and accessories must be dried directly after cleaning.
- Angle rotors, container and hanger made of aluminium are to be lightly greased after drying using acid-free grease, e.g. vaseline.
- The sealing rings of bio-safety systems must be cleaned weekly.
The sealing rings are made of silicone. To guarantee the leak-tightness of the bio-safety systems, the sealing rings must not be handled with talcum powder after cleaning or autoclaving.
Each time before using the bio-safety system, all parts of the bio-safety system must be visually inspected for damage. In addition, the sealing ring(s) of the bio-safety system must be checked to make sure they are in the correct installation position.
Damaged parts of the bio-safety system must be exchanged immediately.
If there are signs of crack formation, brittleness or wear, immediately replace the sealing ring in question. In the case of lids with sealing rings which cannot be replaced, the entire lid must be exchanged.
For the available bio-safety systems, see the chapter "Appendix, Rotors and accessories".
- In order to prevent corrosion as a result of moisture between the rotor and the motor shaft, the rotor should be disassembled and cleaned at least once a month, and the motor shaft should be lightly greased.
- The rotors and accessories have to be checked weekly for wear and corrosion. For swing-out rotors, it is important to check the area of the lifting lugs, for hangers, the grooves and the base should be checked for cracks.

Example: Crack in the groove area:

Rotors and attachments may no longer be utilised upon indication of wear and tear or corrosion.

- Check the firm seating of the rotor on a weekly basis.

29.2.2 Disinfection

- If infectious material should get on the rotors or accessories, they must be appropriately disinfected.
- Ingredients of suitable disinfectants: ethanol, n-propanol, ethyl hexanol, anionic tensides, corrosion inhibitors.
- After using disinfectants, remove disinfectant residue by rinsing with water (only outside of the centrifuge) or wipe off with a damp cloth.
- The rotors and accessories must be dried directly after disinfection.

29.2.3 Removal of radioactive contaminants

- The agent must be specifically labelled as being an agent for the removal of radioactive contaminants.
- Ingredients of suitable agents for removing radioactive contaminants: anionic tensides, non-ionic tensides, polyhydrated ethanol.
- After removing the radioactive contaminants, remove agent residue by rinsing with water (only outside of the centrifuge) or wipe off with a damp cloth.
- The rotors and accessories must be dried directly after removing the radioactive contaminants.

29.2.4 Trunnions

With swing-out rotors the trunnions must be regularly lubricated (Hettich Lubricating Grease No. 4051) in order to ensure consistent swinging out of the hangers.

29.2.5 Rotors and accessories with limited service lives

The use of certain rotors, hangers and accessory parts is limited by time.

These are marked with the maximum permitted number of operating cycles or with an expiration date and the maximum permitted number of operating cycles or just with the expiration date; e.g.:

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For safety reasons, rotors, hangers and accessory parts may no longer be used if either the indicated maximum number of operating cycles or the indicated expiration date has been reached.

29.3 Autoclaving

The following accessories may be autoclaved at 121°C / 250°F (20 min):

- Swing-out rotors
- Angle rotors made of aluminium
- Hangers made of metal
- Lid with bio-seal
- Adapter

Nothing definitive can be said about the degree of sterility.

The lids of the rotors and containers must be removed before autoclaving.

Autoclaving accelerates the ageing process of plastics. In addition, it can cause discolourations in plastics.

The lid of the rotors 1515 and 1515-A may only be autoclaved 10x. Afterwards, it must be exchanged for safety reasons.

After autoclaving, the rotors and the accessories must be visually inspected for damage and any damaged parts must be exchanged immediately.

If there are signs of crack formation, brittleness or wear, immediately replace the sealing ring in question.

In the case of lids with sealing rings which cannot be replaced, the entire lid must be exchanged.

To guarantee the leak-tightness of the bio-safety systems, the sealing rings must not be handled with talcum powder after autoclaving.

29.4 Centrifuge containers

- With leakiness or after the breakage of centrifuging containers broken container parts, glass splinters and leaked centrifugation material are to be completely removed.
- The rubber inserts as well as the plastic sleeves of the rotors are to be replaced after a glass breakage.

Remaining glass splinters cause further glass breakage!

- If this concerns infectious material, a disinfection process is to be executed immediately.

30 Faults

If the fault cannot be eliminated with the help of the fault table, please inform Customer Service.

Please specify the type of centrifuge and the serial number. Both numbers can be found on the name plate of the centrifuge.

Perform a MAINS RESET:

- Switch off the mains switch (switch position "0").
- Wait at least 10 seconds and then switch on the mains switch again (switch position "I").

Message / fault		Cause	Remedy
No display	---	No voltage. Overvoltage protection tripped out.	– Check supply voltage. – Mains switch ON.
TACHO - ERROR	1, 2, 96	Faulty speedometer. Motor, electronics defective.	– Open the cover. – Switch off the mains switch (switch position "0"). – Wait at least 10 seconds. – Turn the rotor vigorously by hand. – Switch on the mains switch again (switch position "I"). The rotor must turn during switch-on.
CONTROL - ERROR	8	Error in lid locking or lid closure.	– Open lid. – Check the loading of the rotor, see chapter "Loading the rotor". – Repeat the centrifugation run.
IMBALANCE	---	The rotor is unevenly loaded.	– Open lid. – Check the loading of the rotor, see chapter "Loading the rotor". – Repeat the centrifugation run.
CONTROL - ERROR	4, 6	Error in lid locking or lid closure.	– Perform a MAINS RESET.
N > MAX	5	Rotation too fast	
N < MIN	13	Rotation too slow	
MAINS INTERRUPT	---	Power failure during the centrifugation run. (The centrifugation run was not finished.)	– Open lid. – Push START / IMPULS button. – Repeat the centrifugation run if necessary.
ROTORCODE	10.1, 10.2	Incorrect rotor coding	– Open lid.
CONTROL-ERROR	21, 22, 25, 27, 29	Error / defect electronics	– Perform a MAINS RESET.
CONTROL-ERROR	23	Error / defect controls	
SER I/O - ERROR	30, 31, 33, 36	Error / defect electronics	
° C * - ERROR	51 - 53, 55	Error / defect electronics	
FU / CCI - ERROR	60 - 64, 67, 68, 82 - 86	Error / defect electronics / motor	
SYNC-ERROR	90	Error / defect electronics	
SENSOR-ERROR	91 - 93	Error / defect unbalance sensor	
KEYBOARD-ERROR	---	Error / defect controls	
NO ROTOR	---	No rotor installed	– Open lid. – Install rotor.
N > ROTOR MAX	---	Speed in the selected program greater than the maximum speed of the rotor. The rotor has been changed. The maximum speed of the installed rotor is higher than that of the rotor that was previously used and this has not yet been detected by rotor recognition.	– Check the set speed. Reduce the set speed – Set a speed up to the maximum speed of the rotor that was previously used. Press the START / IMPULS key to run rotor recognition. See chapter, "Rotor Recognition".

31 Returning Devices

Before returning the device, a transport securing device has to be installed.

If the device or its accessories are returned to Andreas Hettich GmbH & Co. KG, in order to provide protection for people, the environment and materials, it has to be decontaminated and cleaned before being shipped.

We reserve the right to refuse contaminated devices or accessories.

Costs incurred for cleaning and disinfection are to be charged to the customer.

We ask for your understanding in this matter.

32 Disposal

Before disposal, the device must be decontaminated and cleaned to protect people, the environment and property.

When you are disposing of the device, the respective statutory rules must be observed.

Pursuant to guideline 2002/96/EC (WEEE), all devices supplied after August 13, 2005 may not be disposed as part of domestic waste. The device belongs to group 8 (medical devices) and is categorized in the business-to-business field.

The icon of the crossed-out trash can shows that the device may not be disposed as part of domestic waste.

The waste disposal guidelines of the individual EC countries might vary. If necessary, contact your supplier.

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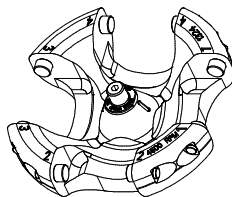
2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)

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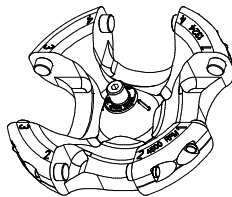
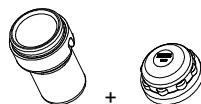
4) can not be centrifugated when plug is attached, Scal. 10µl-300µl, 15ml, 30ml

5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".

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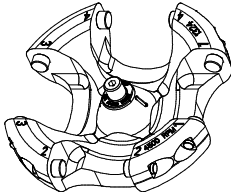
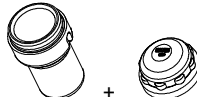
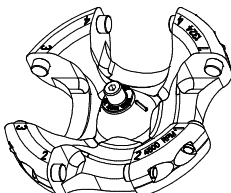
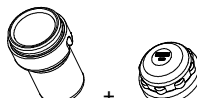
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".

1324	1490 + 1492								
Ausschwingrotor 4-fach / Swing out rotor 4-times  ∠ 90°	 mit Bioabdichtung / with bio-containment ⁵⁾								
	1383					1384	1396	1457	
									
						0513 	0547 		
	Kapazität / capacity	ml	2,6 – 2,9	4,9	4,5 - 5	1,6 -5	4 –7	50	85
Maße / dimensions Ø x L	mm	13 x 65	13 x 90	11 x 92	13 x 75	13 x 100	29 x 115	38 x 106	8 x 66
Anzahl p. Rotor / number p. rotor		20	20	20	20	20	4	4	28
Drehzahl / speed	RPM	4500	4500	4500	4500	4500	4500	4500	4500
RZB / RCF	³⁾	3192	3192	3192	3129	3129	3328	3260	3215
Radius / radius	mm	141	141	141	141	141	147	144	142
 9 (97%)	sec	27							
 9	sec	30							
Temperatur / temperature	°C ¹⁾	- 6							
Probenerwärmung/Sample temp. rise	K ²⁾	10							

1324	1490 + 1492								
Ausschwingrotor 4-fach / Swing out rotor 4-times  ↙ 90°	 mit Bioabdichtung / with bio-containment ⁵⁾								
	1459		4416	4417	6311	6318	1356	0761	
									
			0546	0545			0509	0526	
	 						 ⁷⁾		
Kapazität / capacity	ml	4 – 5,5	7,5 – 8,2	50	30	12	50	15	100
Maße / dimensions Ø x L	mm	15 x 75	15 x 92	29 x 107	26 x 95	17 x 100	29 x 115	17 x 120	44 x 100
Anzahl p. Rotor / number p. rotor		16	16	4	4	4	4	12	4
Drehzahl / speed	RPM	4500	4500	4500	4500	4500	4500	4500	4500
RZB / RCF	³⁾	3215	3215	3283	3056	3328	3328	3328	3192
Radius / radius	mm	142	142	145	135	147	147	147	141
 9 (97%)	sec	27							
 9	sec	30							
Temperatur / temperature	°C ¹⁾	- 6							
Probenerwärmung/Sample temp. rise	K ²⁾	10							

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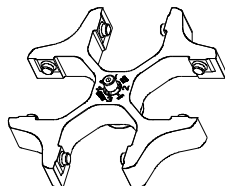
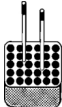
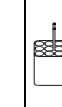
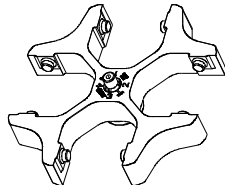
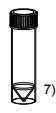
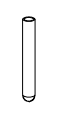
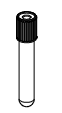
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
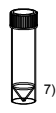
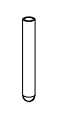
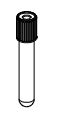
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
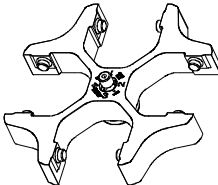
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
 3) Observe the tube manufacturer's instructions.

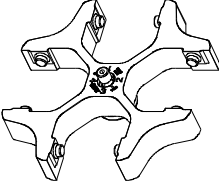
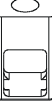
1494	1427 + 1421																
Ausschwingrotor 4-fach / Swing out rotor 4-times  	 + 																
	1357		5229 ⁷⁾		5229		1326		5230		5230		5231		5231 ⁷⁾		
																	
	Rhe- sus	---					0501		0578				0507	0518			
																	
Kapazität / capacity	ml	1	0,4	4,5 – 5		2,7 - 3		6	4	7		4,5 – 5		15		8,5 - 10	
Maße / dimensions $\varnothing \times L$	mm	6 x 45		11 x 92		11x 66		12×82	12×60	12 x 100		11 x 92		17 x 100		16 x 100	
Anzahl p. Rotor / number p. rotor		120		48		48		48		48		48		24		24	
Drehzahl / speed	RPM	5000		5000		5000		5000		5000		5000		5000		5000	
RZB / RCF	³⁾	4081		4053		4053		4053		3941		3941		3941		3941	
Radius / radius	mm	146		145		145		145		141		141		141		141	
 9 (97%)	sec	30															
 9	sec	32															
Temperatur / temperature	$^{\circ}\text{C}$ ¹⁾	- 7															
Probenerwärmung/Sample temp. rise	K ²⁾	+ 15															

1494	1427 + 1421								
Ausschwingrotor 4-fach / Swing out rotor 4-times  	 + 								
	1731	1732			1732		5237	5279	
									
							0500		
									
Kapazität / capacity	ml	25	5	1,6 – 5	2,6 – 2,9	4,9	4 – 7	9	4 – 5,5
Maße / dimensions $\varnothing \times L$	mm	25 x 90	12/13x75	13 x 75	13 x 65	13 x 90	13 x 100	14 x 100	15 x 75
Anzahl p. Rotor / number p. rotor		8	32	32	32	32	32	24	20
Drehzahl / speed	RPM	5000	5000	5000	5000	5000	5000	5000	5000
RZB / RCF	³⁾	3969	4025	4025	4025	4025	4025	3941	4109
Radius / radius	mm	142	144	144	144	144	144	141	147
 9 (97%)	sec	30							
 9	sec	32							
Temperatur / temperature	°C ¹⁾	- 7							
Probenerwärmung/Sample temp. rise	K ²⁾	15							

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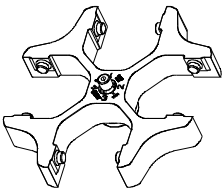
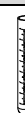
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
3) Observe the tube manufacturer's instructions.
7) not possible to close the lid

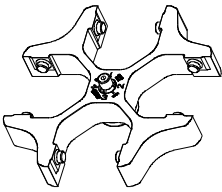
1494		1427 + 1421								
Ausschwingrotor 4-fach / Swing out rotor 4-times   90°		 + 								
		5279	5278	1327	5233	5271		5232	5275 ⁷⁾	
										
					0521	 ⁶⁾		0519	0509	
		Kapazität / capacity	ml	7,5 – 8,2	1,1 – 1,4	3	50	9 - 10	4 - 7	25
Maße / dimensions	Ø x L	mm	15 x 92	8 x 66	10 x 60	34 x 100	16 x 92	16 x 75	24 x 100	17 x 120
Anzahl p. Rotor / number p. rotor		20	48	48	4	20	20	8	4	
Drehzahl / speed	RPM	5000	5000	5000	5000	5000	5000	5000	5000	
RZB / RCF	³⁾	4109	4109	4053	3941	3969	3969	3941	4165	
Radius / radius	mm	147	147	145	141	142	142	141	149	
 9 (97%)	sec	30								
 9	sec	32								
Temperatur / temperature	°C ¹⁾	- 7								
Probenerwärmung/Sample temp. rise	K ²⁾	+ 15								

1494		1427 + 1421				1425			
Ausschwingrotor 4-fach / Swing out rotor 4-times   90°		 + 							
		5276	5277	5272	5273	1432	1433	1434	
									
		0513	2078	0536	---	Rhe-	---	0501	0578
									
Kapazität / capacity	ml	50	1,5	2,0	30	1,5	2,0	1	0,4
Maße / dimensions	Ø x L mm	29 x 115	11 x 38	25 x 110	11 x 38	6 x 45	10 x 60	12 x 82	12 x 100
Anzahl p. Rotor / number p. rotor		4	36	4	48	144	56	48	48
Drehzahl / speed	RPM	5000	5000	5000	5000	5000	5000	5000	5000
RZB / RCF	³⁾	4053	4109	4025	4053	3969	3913	3913	3913
Radius / radius	mm	145	147	144	145	142	140	140	140
 9 (97%)	sec	30							
 9	sec	32							
Temperatur / temperature	°C ¹⁾	- 7							
Probenerwärmung/Sample temp. rise	K ²⁾	15							

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- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
 3) Observe the tube manufacturer's instructions.
 7) not possible to close the lid
 6) Remove the inserts

1494	1425								
Ausschwingrotor 4-fach / Swing out rotor 4-times  ∠ 90°									
	1431				1436	1437	1458	1435	1439
									
	0500	0507	0518		0521	0526		0519	
									
Kapazität / capacity	ml	9	15	7,5 - 8,2	50	100	1,1 – 1,4	25	9 – 10
Maße / dimensions $\varnothing \times L$	mm	14 x 100	17 x 100	15 x 92	34 x 100	44 x 100	8 x 66	24 x 100	16 x 92
Anzahl p. Rotor / number p. rotor		28	28	28	4	4	36	8	16
Drehzahl / speed	RPM	5000	5000	5000	5000	5000	5000	5000	5000
RZB / RCF	³⁾	3913	3913	3913	3913	3801	4025	3913	3913
Radius / radius	mm	140	140	140	140	136	144	140	140
 9 (97%)	sec	30							
 9	sec	32							
Temperatur / temperature	°C ¹⁾	- 10							
Probenerwärmung/Sample temp. rise	K ²⁾	12							

1494	1425								
Ausschwingrotor 4-fach / Swing out rotor 4-times  									
	1438								
									
									
Kapazität / capacity	ml	2,6 – 2,9	2,7 – 3	4,9	4,5 – 5	1,6 – 5	4 – 7	5	4
Maße / dimensions $\varnothing \times L$	mm	13 x 65	11 x 66	13 x 90	11 x 92	13 x 75	13 x 100	12/13x75	12 x 60
Anzahl p. Rotor / number p. rotor		28	28	28	28	28	28	28	28
Drehzahl / speed	RPM	5000							
RZB / RCF	³⁾	3913							
Radius / radius	mm	140							
 9 (97%)	sec	30							
 9	sec	32							
Temperatur / temperature	°C ¹⁾	- 10							
Probenerwärmung/Sample temp. rise	K ²⁾	12							

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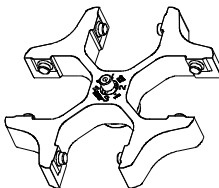
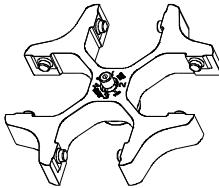
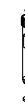
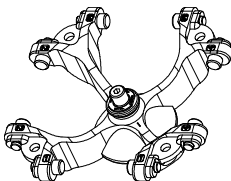
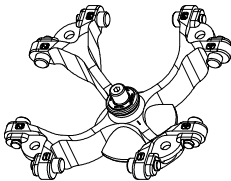
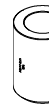
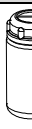
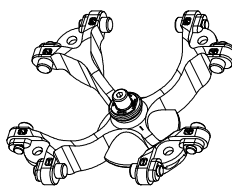
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 4) can not be centrifugated when plug is attached, Scal. 10µl-300µl, 15ml, 30ml
- 5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".

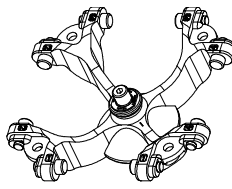
1554	1560 + 1561 / 1565									
Ausschwingrotor 4-fach / Swing out rotor 4-times  ∠ 90° max. Laufzyklen / max. cycles 100 000	 +  /  mit Bioabdichtung / with bio-containment 5) max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 290 g									
	1571		1593	1589			1588			
										
	2078	0536	----	0553	0501	0578	0500	0507	0518	
										
	Kapazität / capacity	ml	1.5	2.0	5	5	6	7	9	15
Maße / dimensions Ø x L	mm	11 x 38		17 x 59	12 x 75	12 x 82	12 x 100	14 x 100	17 x 100	
Anzahl p. Rotor / number p. rotor		56		16	28			20		
Drehzahl / speed	RPM	4500								
RZB / RCF	3)	3328/2332		3328	3215					
Radius / radius	mm	147/103		147	142					
 9 (97%)	sec	28								
 9	sec	31								
Temperatur / temperature	°C 1)	- 8								
Probenerwärmung/Sample temp. rise	K 2)	10								

1554	1560 + 1561 / 1565								
Ausschwingrotor 4-fach / Swing out rotor 4-times  ↙ 90° max. Laufzyklen / max. cycles 100 000	 mit Bioabdichtung / with bio-containment 5) max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 290 g								
	1572	1573	1574	1575	1576	1594		---	
									
	Corex®	0519	0521	0549	0526	---			
							 7)		
Kapazität / capacity	ml	15	25	50	85	100	125	150	200
Maße / dimensions $\varnothing \times L$	mm	18 x 102	24 x 100	34 x 100	38 x 101	44 x 100	51 x 100	51 x 116	56 x 112
Anzahl p. Rotor / number p. rotor		16	4						
Drehzahl / speed	RPM	4500							
RZB / RCF	3)	3260	3056	3124	3260	3124	3328		
Radius / radius	mm	144	135	138	144	138	147		
 9 (97%)	sec	28							
 9	sec	31							
Temperatur / temperature	°C 1)	- 8							
Probenerwärmung/Sample temp. rise	K 2)	10							

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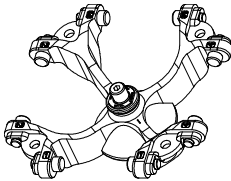
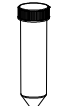
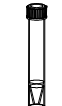
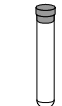
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".
- 7) not possible to close the lid

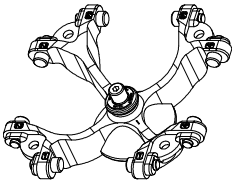
1554	1560 + 1561 / 1565										
Ausschwingrotor 4-fach / Swing out rotor 4-times  90° max. Laufzyklen / max. cycles 100 000	 mit Bioabdichtung / with bio-containment ⁵⁾ max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 290 g										
	1589					1588					
											
											
	Kapazität / capacity	ml	1,1 - 1,4	2,6 - 3,4	4,9	2,7 - 3	4 - 5,5	4 - 5,5	7,5 - 8,5	9 - 10	10
	Maße / dimensions Ø x L	mm	8 x 66	13 x 65	13 x 90	11 x 66	11 x 92	15 x 75	15 x 92	16 x 92	15 x 102
Anzahl p. Rotor / number p. rotor		28					20				
Drehzahl / speed	RPM	4500									
RZB / RCF	³⁾	3215									
Radius / radius	mm	142									
 9 (97%)	sec	28									
 9	sec	31									
Temperatur / temperature	°C ¹⁾	- 8									
Probenerwärmung/Sample temp. rise	K ²⁾	10									

1554	1560 + 1561 / 1565							
Ausschwingrotor 4-fach / Swing out rotor 4-times   max. Laufzyklen / max. cycles 100 000	 mit Bioabdichtung / with bio-containment 5) ----- max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 290 g							
	1589		1588		1591	1581 + E2109	1577	1595
								
	---	---	---	---	---	0509		
								
								7)
Kapazität / capacity	ml	1,6 - 5	4 - 7	4 - 7	8,5 - 10	12	11	15
Maße / dimensions $\varnothing \times L$	mm	13 x 75	13 x 100	16 x 75	16 x 100	17 x 102	16 x 110	17 x 120
Anzahl p. Rotor / number p. rotor		28		20		12	16	8 12
Drehzahl / speed	RPM	4500						
RZB / RCF	3)	3215					3260	3328
Radius / radius	mm	142					144	147
 9 (97%)	sec	28						
 9	sec	31						
Temperatur / temperature	°C 1)	- 8						
Probenerwärmung/Sample temp. rise	K 2)	10						

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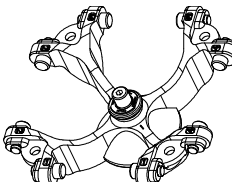
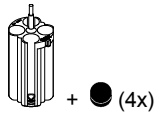
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 5) In conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".
- 7) not possible to close the lid

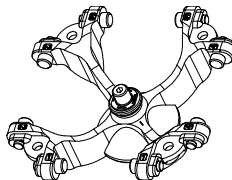
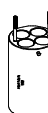
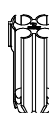
1554	1560 + 1561 / 1565								
Ausschwingrotor 4-fach / Swing out rotor 4-times  $\angle 90^\circ$ max. Laufzyklen / max. cycles 100 000	 +  /  mit Bioabdichtung / with bio-containment 5) ----- max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 290 g								
	1578	1579	1581	1582	1583	1584	1585		
									
	----	0513	----	----	----		0545		
									
Kapazität / capacity	ml	30	50	12	25	30	50	10	30
Maße / dimensions $\varnothing \times L$	mm	25 x 110	30 x 115	17 x 100	25 x 90	25 x 110	29 x 115	16 x 80	26 x 95
Anzahl p. Rotor / number p. rotor		4		16		4		20	4
Drehzahl / speed	RPM	4500							
RZB / RCF	3)	3328		3260		3328		3215	3260
Radius / radius	mm	147		144		147		142	144
 9 (97%)	sec	28							
 9	sec	31							
Temperatur / temperature	°C 1)	- 8							
Probenerwärmung/Sample temp. rise	K 2)	10							

1554	1560 + 1561 / 1565						
Ausschwingrotor 4-fach / Swing out rotor 4-times  $\angle 90^\circ$ max. Laufzyklen / max. cycles 100 000	 mit Bioabdichtung / with bio-containment 5) -----						
	max. Laufzyklen / max. cycles: 80000						
	max. Beladung / max. load: 290 g						
	1586	1575	1587				
							
Kapazität / capacity	ml	50	85	94	30		
Maße / dimensions $\varnothing \times L$	mm	29 x 107	38 x 106	38 x 102	44 x 105		
Anzahl p. Rotor / number p. rotor		4					
Drehzahl / speed	RPM	4500					
RZB / RCF	³⁾	3260		3192			
Radius / radius	mm	144		141			
 9 (97%)	sec	28					
 9	sec	31					
Temperatur / temperature	$^{\circ}\text{C}^{1)}$	- 8					
Probenerwärmung/Sample temp. rise	K ²⁾	10					

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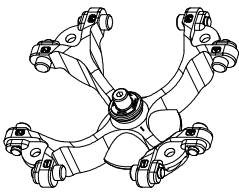
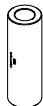
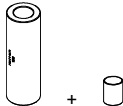
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 4) can not be centrifugated when plug is attached, Scal. 10µl-300µl, 15ml, 30ml
- 5) In conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".

1554				1559														
Ausschwingrotor 4-fach / Swing out rotor 4-times  $\angle 90^\circ$ max. Laufzyklen / max. cycles 100 000				 max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 200 g														
				1486			1482A			1482A + 4 x 0715			1486			1482A		
																		
				0553	0501	0578	0500	0507	0518									
																		
Kapazität / capacity	ml	5	6	7	9	15	4 - 7	4 - 5,5	2,6 - 3,4	2,7 - 3	4,5-5	4,9	9-10	10				
Maße / dimensions $\varnothing \times L$	mm	12 x 75	12 x 82	12 x 100	14 x 100	17 x 100	16 x 75	15 x 75	13 x 65	11 x 66	11 x 92	13 x 90	16 x 92	15 x 102				
Anzahl p. Rotor / number p. rotor		20			16			16		20		20		16				
Drehzahl / speed	RPM	4500																
RZB / RCF	³⁾	3215			3260			3215		3215			3260					
Radius / radius	mm	142			144			142		142			144					
 9 (97%)	sec	28																
 9	sec	31																
Temperatur / temperature	°C ¹⁾	- 8																
Probenerwärmung / Sample temp. rise	K ²⁾	9																

1554		1559																				
Ausschwingrotor 4-fach / Swing out rotor 4-times  $\angle 90^\circ$ max. Laufzyklen / max. cycles 100 000	 max. Laufzyklen / max. cycles: 80000 max. Beladung / max. load: 200 g																					
	1486		1482A		1488		1487		1483A		1484		1482A		1484							
																						
									0509		0513		---		---							
																						
	Kapazität / capacity		ml		1,6 - 5		4 - 7		8,5-10		8		12		15		50		12		50	
	Maße / dimensions $\varnothing \times L$		mm		13 x 75		13 x 100		16x100		16 x 125		17 x 102		17 x 120		29 x 115		17 x 100		29 x 115	
Anzahl p. Rotor / number p. rotor		20		16		12		16		4		16		4								
Drehzahl / speed		RPM		4500																		
RZB / RCF		³⁾		3215		3260		3328		3147		3351		3305		3260		3305				
Radius / radius		mm		142		144		147		139		148		146		144		146				
 9 (97%)		sec		28																		
 9		sec		31																		
Temperatur / temperature		°C ¹⁾		- 8																		
Probenerwärmung / Sample temp. rise		K ²⁾		9																		

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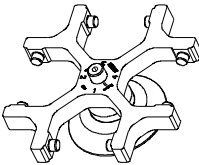
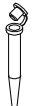
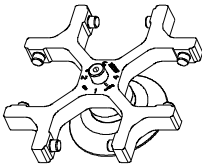
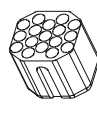
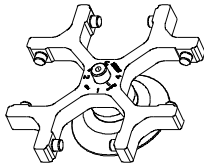
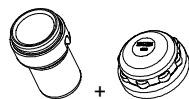
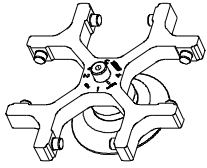
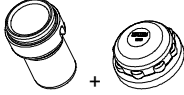
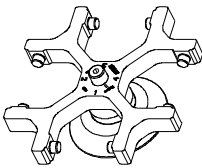
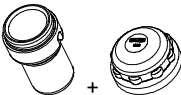
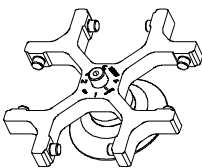
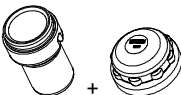
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".
- 6) Remove the inserts

1624	1481 + 1492							
Ausschwingrotor 4-fach / Swing out rotor 4-times   90°	 mit Bioabdichtung / with bio-containment ⁵⁾							
	1383							
								
	0501	0578						
Kapazität / capacity ml	6	7	4,9	4,5 - 5,0	2,7 - 3,0	2,6 - 2,9	1,6 - 5,0	5
Maße / dimensions Ø x L mm	12 x 82	12 x 100	13 x 90	11 x 92	11 x 66	13 x 65	13 x 75	12 x 75
Anzahl p. Rotor / number p. rotor	20	20	20	20	20	20	20	20
Drehzahl / speed RPM	4000	4000	4000	4000	4000	4000	4000	4000
RZB / RCF ³⁾	2558	2558	2558	2558	2558	2558	2558	2558
Radius / radius mm	143	143	143	143	143	143	143	143
 9 (97%) sec	20							
 9 sec	25							
Temperatur / temperature °C ¹⁾	- 15							
Probenerwärmung/Sample temp. rise K ²⁾	7							

1624	1481 + 1492							
Ausschwingrotor 4-fach / Swing out rotor 4-times   90°	 mit Bioabdichtung / with bio-containment ⁵⁾							
	1383	1384	1396	1457	1343		1363	1365
								
		0513	0547					
Kapazität / capacity ml	4 - 7,0	50	85	1,1 - 1,4	3	4	25	30
Maße / dimensions Ø x L mm	13 x 100	29 x 115	38 x 106	8 x 66	10 x 60	10 x 88	25 x 90	25 x 110
Anzahl p. Rotor / number p. rotor	20	4	4	28	36	36	4	4
Drehzahl / speed RPM	4000	4000	4000	4000	4000	4000	4000	4000
RZB / RCF ³⁾	2558	2665	2612	2576	2630	2630	2343	2665
Radius / radius mm	143	149	146	144	147	147	131	149
 9 (97%) sec	20							
 9 sec	25							
Temperatur / temperature °C ¹⁾	- 15							
Probenerwärmung/Sample temp. rise K ²⁾	7							

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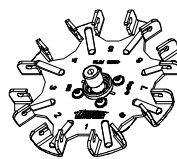
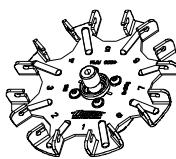
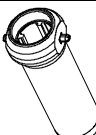
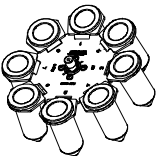
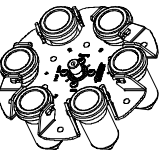
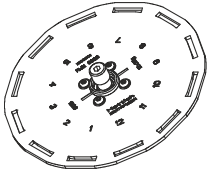
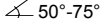
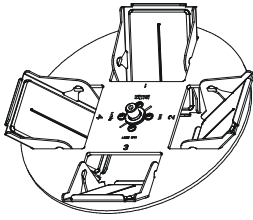
CP Kulturplatte /
Culture plate

DWP Deep Well Platte /
Deep well plate

MS Micronic System /
Micronic system

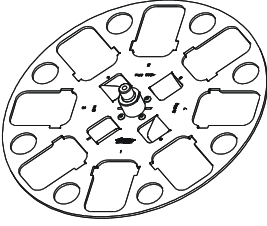
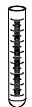
1628	1127 (50°)		1645	4345	
Ausschwingrotor 12-fach / Swing out rotor 12-times  			Ausschwingrotor 4-fach / Swing out rotor 4-times  		
				MTP	CP
					
Kapazität / capacity	ml	2,7 – 3	2,6 – 2,9	Kapazität / capacity	ml
Maße / dimensions Ø x L	mm	11 x 66	13 x 65	Maße / dimensions TxBxH / DxVxH	mm
Anzahl p. Rotor / number p. rotor		12		Anzahl p. Rotor / number p. rotor	4
Drehzahl / speed	RPM	5000		Drehzahl / speed	RPM
RZB / RCF		3466		RZB / RCF	
Radius / radius	mm	124		Radius / radius	mm
 9 (97%)	sec	16		 9 (97%)	sec
 9	sec	16		 9	sec
Temperatur / temperature	°C ¹⁾	- 15		Temperatur / temperature	°C ¹⁾
Probenerwärmung/Sample temp. rise	K ²⁾	14		Probenerwärmung/Sample temp. rise	K ²⁾

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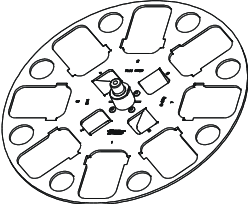
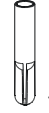
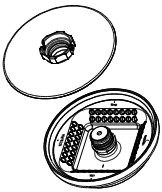
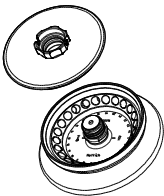
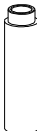
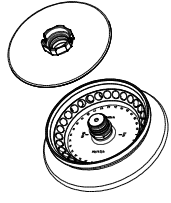
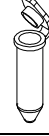
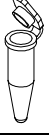
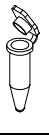
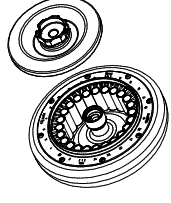
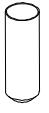
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.
- 5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".
- 8) recommended for high-speed centrifugation
- 9) Load only each second position of the rotor
- 14) With running time longer than 20 min., the sample temp. rise up will be > 20°K (only with cooling centrifuges)

1553								
Winkelrotor 30-fach / Angle rotor 30-times  45° mit Bioabdichtung / with bio-containment ⁵⁾		8) 2031	2023	2024	0788			
								
		0536	2078				---	
								9) 
Kapazität / capacity	ml	2,0	1,5	0,8	0,5	0,4	0,2	0,5
Maße / dimensions	Ø x L mm	11 x 38	11 x 38	8 x 45	8 x 30	6 x 45	6 x 18	10,7 x 36
Anzahl p. Rotor / number p. rotor		30	30	30	30	30	30	15
Drehzahl / speed	RPM	14150	14150	14150	14150	14150	14150	14150
RZB / RCF	³⁾	21713	21713	21713	21713	21713	21713	20818
Radius / radius	mm	97	97	97	97	97	97	93
 9 (97%)	sec	35						
 9	sec	32						
Temperatur / temperature	°C ¹⁾	- 1						
Probenerwärmung/Sample temp. rise	K ¹⁵⁾	20						

1555								
Ausschwingrotor 24-fach / Swing out rotor 24-times  90° mit Bioabdichtung / with bio-containment ⁵⁾		2024	2023	8) 2031				
					---			
		---	---	2078	0536			
								
Kapazität / capacity	ml	0,2	0,4	0,5	0,8	1,5	2,0	
Maße / dimensions	Ø x L mm	6 x 18	6 x 45	8 x 30	8 x 45	11 x 38	11 x 38	
Anzahl p. Rotor / number p. rotor		24						
Drehzahl / speed	RPM	13000						
RZB / RCF	³⁾	18327						
Radius / radius	mm	97						
 9 (97%)	sec	36						
 9	sec	31						
Temperatur / temperature	°C ¹⁾	3						
Probenerwärmung/Sample temp. rise	K ¹⁰⁾	20						

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1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)

3) Observe the tube manufacturer's instructions.

5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio

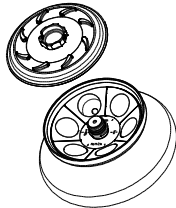
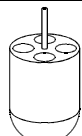
safety systems in chapters "Notes on safety" and "Maintenance and servicing".

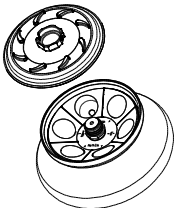
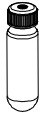
8) recommended for high-speed centrifugation

9) Load only each second position of the rotor

10) With running time longer than 10 min., the sample temp. rise up will be > 20°K (only with cooling centrifuges)

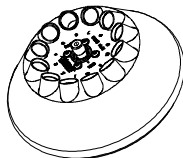
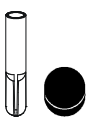
15) With running time longer than 15 min., the sample temp. rise up will be > 20°K (only with cooling centrifuges)

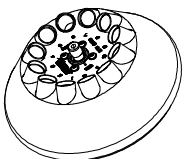
1556											
Winkelrotor 6-fach / Angle rotor 6-times  35° mit Bioabdichtung / with bio-containment ⁵⁾	1449		1477	1478							
											
	2078	0536	----	----	0507	0518					
											
	Kapazität / capacity	ml	1,5	2,0	0,5	10	15	7,5 – 8,2	9 - 10	10	8,5 - 10
	Maße / dimensions Ø x L	mm	11 x 38	10,7 x 36	16 x 80	17 x 100	15 x 92	16 x 92	15 x 102	16 x 100	
	Anzahl p. Rotor / number p. rotor		24	24	12	6					
	Drehzahl / speed	RPM	9000								
	RZB / RCF ³⁾		10324	10324	10414	10052					
	Radius / radius	mm	114	114	115	111					
 9 (97%)	sec	36									
 9	sec	32									
Temperatur / temperature	°C ¹⁾	0									
Probenerwärmung/Sample temp. rise	K ²⁾	20									

1556										
Winkelrotor 6-fach / Angle rotor 6-times  35° mit Bioabdichtung / with bio-containment ⁵⁾			1466	1454	1447	1446			1463	
										
	0547	0539 / 538	0509	0513	0545	0546	0549	0548	0521	
										
Kapazität / capacity	ml	85	94	15	50	30	50	85	50	50
Maße / dimensions Ø x L	mm	38 x 106	38 x 102	17 x 120	29 x 115	26 x 95	29 x 107	38 x 101	35 x 105	34 x 100
Anzahl p. Rotor / number p. rotor		6								
Drehzahl / speed	RPM	9000								
RZB / RCF ³⁾		10595	10052		9690	10142	10595	10414		
Radius / radius	mm	117	111		107	112	117	115		
 9 (97%)	sec	36								
 9	sec	32								
Temperatur / temperature	°C ¹⁾	0								
Probenerwärmung/Sample temp. rise	K ²⁾	20								

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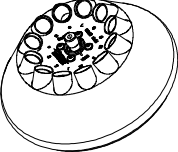
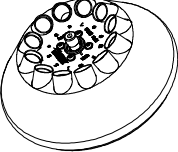
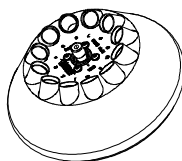
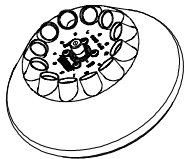
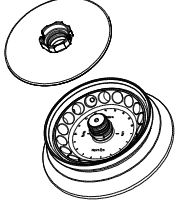
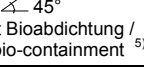
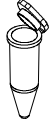
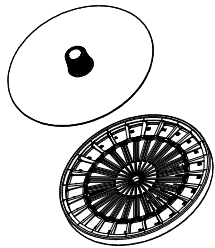
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
3) Observe the tube manufacturer's instructions.

1615									
Winkelrotor 12-fach / Angle rotor 12-times   35°									
		SK 19/85-4							
									
Kapazität / capacity	ml	4,5 - 5	4,9	7,5 – 8,2	9 – 10	10	4 – 5,5	1,6 - 5	4 - 7
Maße / dimensions	Ø x L mm	11 x 92	13 x 90	15 x 92	16 x 92	15 x 102	15 x 75	13 x 75	16 x 75
Anzahl p. Rotor / number p. rotor		12	12	12	12	12	12	12	12
Drehzahl / speed	RPM	12000	12000	12000	12000	12000	12000	12000	12000
RZB / RCF	³⁾	16582	16582	16582	16582	16582	14650	14650	14650
Radius / radius	mm	103	103	103	103	103	91	91	91
 9 (97%)	sec					40			
 9	sec					40			
 0	sec					840			
Temperatur / temperature	°C ¹⁾					- 2			
Probenerwärmung/Sample temp. rise	K ²⁾					14			

1615											
Winkelrotor 12-fach / Angle rotor 12-times  ∠ 35°						6305		1063-6			
						SK 1/89					
			0507	0518	0509	---			2078	0536	
											
Kapazität / capacity	ml	4 - 7	15	15	0,8	4	0,5	1,5	2,0		
Maße / dimensions	Ø x L mm	13 x 100	17 x 100	17 x 120	8 x 45	10 x 88	10,7 x 36	11 x 38			
Anzahl p. Rotor / number p. rotor		12	12	6	12	12	12				
Drehzahl / speed	RPM	12000									
RZB / RCF	³⁾	16582	16582	16582	10303	14006	11108	10947			
Radius / radius	mm	103	103	103	64	87	69	68			
 9 (97%)	sec	40									
 9	sec	40									
Temperatur / temperature	°C ¹⁾	- 2									
Probenerwärmung/Sample temp. rise	K ²⁾	14									

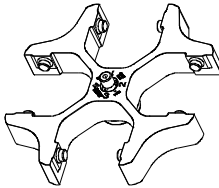
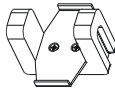
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
- 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
- 3) Observe the tube manufacturer's instructions.

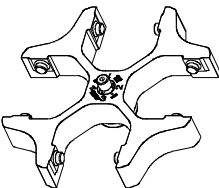
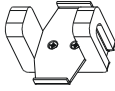
1627										
Winkelrotor 18-fach / Angle rotor 18-times  										
										
Kapazität / capacity	ml	5								
Maße / dimensions Ø x L	mm	17 x 59								
Anzahl p. Rotor / number p. rotor		18								
Drehzahl / speed	RPM	14150								
RZB / RCF	³⁾	22161								
Radius / radius	mm	99								
 9 (97%)	sec	35								
 9	sec	32								
Temperatur / temperature	°C ¹⁾	2								
Probenerwärmung/Sample temp. rise	K ¹⁰⁾	20								

1650-A										
Hämatokritrotor 24-fach / Haematocrite rotor 24-times 										
										
Kapazität / capacity	ml									
Maße / dimensions Ø x L	mm	1,4 x 75								
Anzahl p. Rotor / number p. rotor		24								
Drehzahl / speed	RPM	14150								
RZB / RCF	³⁾	19027								
Radius / radius	mm	85								
 9 (97%)	sec	35								
 9	sec	32								
Temperatur / temperature	°C ¹⁾	- 3								
Probenerwärmung/Sample temp. rise	K ¹⁰⁾	20								

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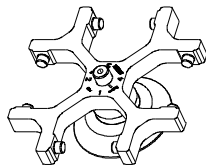
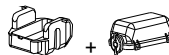
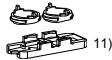
- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
3) Observe the tube manufacturer's instructions.
5) in conformity with DIN EN 61010, part 2 – 020. Observe the notes for bio safety systems in chapters "Notes on safety" and "Maintenance and servicing".
10) With running time longer than 10 min., the sample temp. rise up will be > 20°K (only with cooling centrifuges)

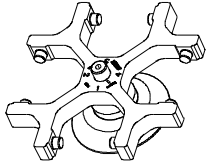
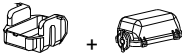
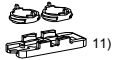
1494	1452								
Ausschwingrotor 4-fach / Swing out rotor 4-times  	1662						1670		
							 11)		
	1663	1664	1665	1666	1667	1668	1663	1664	
									
	Kapazität / capacity	ml	1	2	4	8	3 x 2	4 x 1	1
Maße / dimensions $\varnothing$ / A	mm ²	6,2 / 30	8,7 / 60	12,4 / 120	17,5 / 240	8,7 / 60	6,2 / 30	6,2 / 30	8,7 / 60
Anzahl p. Rotor / number p. rotor		4	4	4	4	4	4	4	4
Filterkarten / filter cards		1675	1675	1675	1676	1677	1678	1692	1692
Drehzahl / speed	RPM	5000							
RZB / RCF	3)	2879							
Radius / radius	mm	103							
 9 (97%)	sec	30							
 9	sec	32							
Temperatur / temperature	°C 1)	- 10							
Probenerwärmung/Sample temp. rise	K 2)	12							

1494	1452							
Ausschwingrotor 4-fach / Swing out rotor 4-times  $\angle 90^\circ$								
	1670				1285	1470		
	 11)							
	1665	1666	1667	1668		1471	1475	
								
Kapazität / capacity	ml	4	8	3 x 2	4 x 1	Objektträger/ object slide	1 x 8	2 x 8
Maße / dimensions $\varnothing$ / A	mm ²	12,4 / 120	17,5 / 240	8,7 / 60	6,2 / 30	26 / 76	17,5 / 240	17,5 / 240
Anzahl p. Rotor / number p. rotor		4	4	4	4	24	4	4
Filterkarten / filter cards		1692	1691	1694	1693	---	---	---
Drehzahl / speed	RPM	5000	5000	5000	5000	5000	5000	5000
RZB / RCF	³⁾	2879	2879	2879	2879	2767	2739	2739
Radius / radius	mm	103	103	103	103	99	98	98
 9 (97%)	sec	30						
 9	sec	32						
Temperatur / temperature	°C ¹⁾	- 10						
Probenerwärmung/Sample temp. rise	K ²⁾	12						

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- 1) Lowest possible temperature during maximum speed, 1 h running time and 20°C ambient temperature (only with cooling centrifuges)
 2) Sample temp. rise during maximum speed and 1 hour running time (only with centrifuges without cooling)
 3) Observe the tube manufacturer's instructions.
 11) Object slide will not stand RCF values exceeding 1100

1624	1660 + 1661								
Ausschwingrotor 4-fach / Swing out rotor 4-times  ∠ 90°									
	1662						1670		
							 11)		
	1663	1664	1665	1666	1667	1668	1663	1664	
									
Kapazität / capacity	ml	1	2	4	8	3 x 2	4 x 1	1	2
Maße / dimensions Ø / A	mm²	6,2 / 30	8,7 / 60	12,4 / 120	17,5 / 240	8,7 / 60	6,2 / 30	6,2 / 30	8,7 / 60
Anzahl p. Rotor / number p. rotor		4	4	4	4	4	4	4	4
Filterkarten / filter cards		1675	1675	1675	1676	1677	1678	1692	1692
Drehzahl / speed	RPM	4000							
RZB / RCF	³⁾	1646							
Radius / radius	mm	92							
 9 (97%)	sec	20							
 9	sec	25							
Temperatur / temperature	°C ¹⁾	- 16							
Probenerwärmung/Sample temp. rise	K ²⁾	3							

1624	1660 + 1661				1660	1680			
Ausschwingrotor 4-fach / Swing out rotor 4-times  $\angle 90^\circ$									
	1670				1285	1662			
	 ¹¹⁾				 nur ohne Deckel 1661 / only without lid 1661				
	1665	1666	1667	1668		1671	1672	1673	
									
Kapazität / capacity	ml	4	8	3 x 2	4 x 1	Objekträger Object slide	[1] 0,5	[1] 0,5	[1] 0,5
Maße / dimensions $\varnothing$ / A	mm ²	12,4 / 120	17,5 / 240	8,7 / 60	6,2 / 30	26 / 76	6,2 / 30	8,7 / 60	12,4 / 120
Anzahl p. Rotor / number p. rotor		4	4	4	4	24	4	4	4
Filterkarten / filter cards		1692	1691	1694	1693	---	[1] 1696	[1] 1697	[1] 1698
Drehzahl / speed	RPM	4000	4000	4000	4000	4000	4000	4000	4000
RZB / RCF	³⁾	1646	1646	1646	1646	1520	1467	1467	1467
Radius / radius	mm	92	92	92	92	88	82	82	82
 9 (97%)	sec	20							
 9	sec	25							
 0	sec	390							
Temperatur / temperature	°C ¹⁾	- 16							
Probenerwärmung/Sample temp. rise	K ²⁾	3							

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All rights reserved. No part of this publication may be reproduced without the prior written permission of the copyright owner.